 Kartell

The image consists of five separate geometric diagrams arranged vertically. From top to bottom: 1) A simple rectangle. 2) A semi-oval shape with a straight line segment tangent to its right side at the top. 3) A triangle with a horizontal line segment passing through its base. 4) A zigzag line consisting of two parallel diagonal line segments. 5) Another semi-oval shape with a straight line segment tangent to its right side at the bottom.

A large, vertical graphic element consisting of the words "The Art of" stacked vertically. The letters are bold and stylized, with the "T" and "A" in "The" being white with red outlines, and the "R" and "t" in "Art" being red with white outlines. The background is a solid blue color.

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Kartell, more than 60 years as leader in industrial design

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KARTELL® LABWARE DIVISION

Kartell, founded in 1949, is one of the first companies in Europe to use plastic materials for the production of design furniture, furnishing accessories and home accessories.

The guideline that has characterized Kartell's over sixty years of activity is the idea of combining the technological and industrial aspect, necessary to respond to specific economic and social needs, to the cultural and aesthetic one. In this sense, Kartell is one of the leading interpreters of industrial design in the panorama of Italian and international companies.

At the end of the 1950s, the evolution in technological-design experimentation leads, with great intuition, to the creation of the Labware Division for the supply of technical articles for laboratories; a division able to make the best use of materials such as polypropylene, polystyrene, polyethylene and to offer plastic as a valid alternative to glass due to its lightness, strength, economy and safety.

In 1958, thanks to the experience of the Housewares Division, it has already gained a good knowledge of industrial production and is open to new development opportunities. Giulio Castelli, a graduate in chemical engineering, has the technical mastery to evaluate these new opportunities wisely.

The chemical-physical specifications, the characteristics of unbreakability and the search for materials, give the new plastics great advantages over glass and make them interesting also for the production of laboratory items such as beakers, cylinders, funnels and test tubes. From the first experiences Giulio Castelli entrusts the design of the Labware objects to his designers, Gino Colombini first and Centrokappa after: a choice that is one of the keys to the success of the Division.

As already done with housewares, Gino Colombini tackles the design of technical articles by tracing innovative formal solutions, with small or large changes to the traditional form of the object that improve its functionality. Proof of this, among others, are the long-stemmed funnels of 1962, with small ribs on the inside to prevent the filter paper from sticking to the surface or the test tube holder from 1959, for which Colombini identifies a new modular solution with "X" comb elements that replace the traditional support with overlapping planes with large holes. Or again, from Centrokappa from 1979, the First EYE first aid in which the support also has a signaling function.

For the realization of its Labware articles, Kartell uses over forty different plastic materials: the Division is a technological gym where the company conducts research on materials and identifies the most suitable technologies for each individual product.

The choice of plastics to be used depends on the different applications of the articles; for this reason in the production process there are important elements such as concentrations and times of contact with chemical agents, aging, thermal stress, aggressive detergents, exposure to U.V. radiation, sterilization and other environmental effects.

Having always favored quality, the use of ever more advanced technologies, the choice of the range of items in the catalog, makes the Kartell image and name recognized worldwide.

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COMPANY QUALITY CERTIFICATION: ISO 9001:2015

In 1996, Kartell decided to certify its Corporate Quality Management System in compliance with UNI EN ISO 9001:1994 standards.

In 2005, the company aligned its Quality Management Systems with the standard UNI EN ISO 9001:2000.

In 2008, the company renewed its ISO 9001:2000 certification.

In 2010 Kartell updated its certification standard and is now UNI EN ISO 9001:2008 certified. A guarantor for this certification process is the I.I.P. (Italian Institute of Plastics), which is itself accredited by SINCERT and CISQ, the Italian federation of accreditation bodies for Quality Management Systems.

In 2017, Kartell updated its certification standard to UNI EN ISO 9001:2015.

CISQ is part of IQNET (International Certification Network), a supranational body which guarantees mutual recognition of the ISO standard in countries worldwide. The attainment and maintenance of this certification, made possible by the commitment and perseverance of all company offices involved, testifies to the continued research into ever higher levels of quality in company management systems. A copy of the Quality Certification can be downloaded from the corporate website www.kartelllabware.com

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CERTIFICATE

CISQ/IIP S.r.l.
as an IQNet Partner hereby states that the organization:

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VIA DELLE INDUSTRIE, 1 - 20082 NOVIGLIO (MI)

for the following scope:

Design, industrialization and production of furnishing accessories. Design and production of laboratory technical items.

has implemented and maintains a

Quality Management System

which fulfills the requirements of the following standard

ISO 9001:2015

Issued on: 2017/07/28

First issued on: 2005/03/14
for the validity date, please refer to the original certificate* issued by CISQ/IIP

Registration Number:
IT-16397 IIP 694

Michael Drechsel
President of IQNET

Ing. Claudio Proveti
President of CISQ

AENOR Spain AFNOR Certification France Vlaamse Belgische APCER Portugal CCC Cyprus
CGC Italy CGC Italia CONCAVE CO CDS Czech Republic Cetis Chile DQS Holding GmbH Germany
FCAV Argentina IQA Japan KFO Korea MIRTec Greece MSZT Hungary NemoK AS Norway NSAI Ireland PCBC Poland
Quality Austria RIA Russia SIGE Mexico SH Israel SIQ Slovenia SIRIM QAS International Malaysia
SQS Switzerland SRAC Romania TEST St Petersburg Russia TSE Turkey YUQS Serbia
TUV Rheinland Turkey TÜV SÜD Switzerland UNEVOC United Arab Emirates UQAS UAE Inc.

* This attestation is directly linked to the IQNet Partner's original certificate and shall not be used as a stand-alone document

** The list of IQNet partners is valid at the time of issue of this certificate. Updated information is available under www.iqnet-certification.com

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ACCREDIA
SGD N° 200A
SGC N° 0008
SGR N° 014P
LAB N° 0007
Memorandum of Understanding
Accreditation of Laboratory
Signature of EA, IAF and ILAC Mutual Recognition Agreement

IAP SRL - via Veltella 2 - 20900 Monza (MB)
www.iap.it info@iap.it

CISQ is a member of
IQNet, the association of the world's leading
certification bodies for management systems.
IQNet is composed of more than 30 bodies and covers
over 150 countries of Latin America.

CISQ is the Italian Federation of Management
Systems Certification Bodies.

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CERTIFICATIONS: ISO 14001:2015

RECYCLABILITY

Recyclability, sustainability, eco-compatibility - in a nutshell, respect for the environment - these are the themes which the world community takes very much to heart. It may seem absurd to talk about environment in terms of productions having plastic materials as their common denominator. Nevertheless, contrary to common belief, Kartell products are perfectly recyclable. To simplify the recycling process, in fact, the various components of Kartell products can be easily separated out and traced back to their single material elements. The plastic parts of each product also bear clear identification markings in accordance with the DIN 6120 specification, part 2, just so that they can be correctly identified and to facilitate recycling.

DIN 6120 PART 2 - RECYCLING SYMBOLS

The code within the logo identifies the compound in accordance with the following table:

01	02	03	04	05	06	07
PET	PE-HD	PVC	PE-LD	PP	PS	0

ISO 14001 CERTIFICATION: 2004

In 2011 Kartell achieved UNI EN ISO 14001:2004 certification for its support of an effective Environmental Management System, an international standard recognised throughout the world and developed about 10 years ago which defines development and implementation parameters in corporate processes in order to achieve an effective environmental management system.

WHAT IS ISO 14001: 2004?

This certification attests that the organization certified has implemented a management system capable of controlling environmental impact in its own business and systematically endeavours to improve it in a sustainable, effective and consistent manner. ISO 14001 certification is not obligatory but is the result of the voluntary choice of the Company which decides to define, implement, maintain and improve its own environmental management system. The production of high quality, safe and reliable design products with low environmental impact and which are recyclable. This is the challenge that Kartell is called upon to come to grips with as protagonist in the coming years.

In 2017, Kartell updated the certification standard to UNI EN ISO 14001:2015. A copy of the Quality Certification is available for download on www.kartelllabware.com

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THE INTERNATIONAL CERTIFICATION NETWORK

CERTIFICATE

CISQ/IIP S.r.l.

as an IQNet Partner hereby states that the organization:

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VIA DELLE INDUSTRIE, 1 - 20082 NOVIGLIO (MI)

for the following scope:

Design, production management of furnishing and design accessories through the transformation of plastics, metal and wood materials.

Design and production management of technical laboratory items through the transformation of plastic materials.

has implemented and maintains an

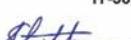
Environmental Management System

which fulfills the requirements of the following standard

ISO 14001:2015

Issued on: 2018/01/12
First issued on: 2012/01/13
for the validity date, please refer to the original certificate issued by CISQ/IIP

Registration Number:
IT-58594 IIP 170

 
Alex Stochitoiu
President of IQNET

IQNet Partners**:

AFOR Spain AFNOR Certification France APCER Portugal CCC Cyprus CISQ Italy

CCS China CQC China CCC Registration Bureau Cen-Croatia DQS Holding GmbH Germany FCAS Brazil

FONDRONOMA France FQS Czech Republic GRCO Colombia GRS Certification Foundation INTECO Costa Rica

Quality ROMA Austria KPF Korea MIRTIC Greece MSCT Hungary Nemetsz AS Norway NSAI Ireland PCBC Poland

SQS Switzerland SIRIM Malaysia SII Israel SQS Slovakia SIRIM QAS International Malaysia

SQS Switzerland SIRIM Malaysia SII Israel SQS Slovakia SIRIM QAS International Malaysia

SGS Switzerland SWAC Romania TEST Testing Services Russia TUV Rheinland Germany TÜV Nord Germany

TÜV Nord Germany TÜV Nord Germany TÜV Nord Germany TÜV Nord Germany TÜV Nord Germany

** This annotation is directly linked to the IQNet Partner's original certificate and shall not be used as a stand-alone document

** The list of IQNet partners is valid at the time of issue of this certificate. Updated information is available under www.iqnet-certification.com

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Design, production management of furnishing and design accessories through the transformation of plastics, metal and wood materials.
Design and production management of technical laboratory items through the transformation of plastic materials.

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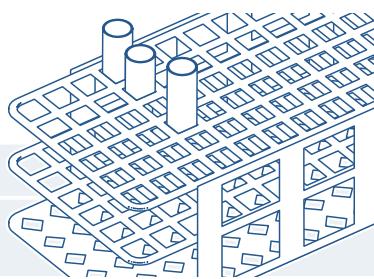
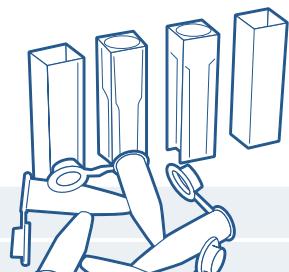
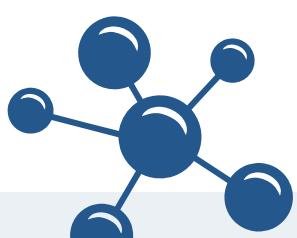
ALPHABETICAL INDEX

#

4 Way connector	17
25 Slides staining trough	59
120° "Y" tapered connectors	17
A	
Adhesive labels	42
Adjustable volume dispenser bottles	43
Analytical funnels	67
Aspirator bottles, heavy duty	45
B	
Balance boats	128
Base plate for test tube racks	81
Beakers	95
Bed pan	87
Bench pipette holder	72-77
Blue graduated beakers, low form - class b	50
Blue graduated measuring cylinders tall form - class b	48-49
Boehringer es 300®, es 600® sample cup	103
Bottle-carriers	43
Bottles	95
Bucket with spout	31
Burette clamps	71
C	
Cap for cuvettes	101
Capsules	95
Centrifichem®, technicon®, beckman® sample cup	104
Centrifuge tubes, cylindrical	80
Certifications: iso 14001:2015	5
Clamps for glass joints	71
Cobas mira® multi-cuvettes	101
Colorimetric 8 cell tray	61
Colorimetric tray	61
Company quality certification: iso 9001:2015	4
Conical 1 and 2 hole stoppers	23
Conical centrifuge tubes	80
Conical centrifuge tubes, graduated	80
Conical erlenmeyer flasks	53
Conical measures, graduated	51
Conical stoppers	23-24- 52
Container for urine	87
Containers for urine and biological specimens	109
Coulter® cell counter cup	103
Crystallisation vessels	95
Cuvette for olli®	104
Cuvettes for amelung®	105
Cuvettes for fibrintimer®	105
Cuvettes holder	101
Cylindrical jars with ribbed cap	42-60
Cylindrical test tubes	105
Deep tray, low form, stackable	30
Deep tray, low form with handles and draining holes	31
D	
Deep trays	30
Desiccator plates	64
Disconnectors - tapered	18
Disks for buchner funnels	66
Dispensing bottle	40
Disposable	106
Disposable dropping pipettes - bellow type	125

Disposable petri dishes

126	Hand operated vacuum pump
101	Heavy-walled carboy bottles, narrow neck
106	Heavy-walled carboy bottles, wide neck
132	Hoses
130	Hoses connectors
131	Hoses k30
131	Hoses k70
131	Hoses selector
25	Hot grip
87	I
70	Ice bucket
86	Imhoff cone stand
41	Industrial scoops
25	Industrial tanks
18	Inoculating loops and needles
22	Input pot
127	Input trays
124	Integral eye wash bottle m45
119	Integral wash bottles
110	Invalid cup
98	Isoflex laboratory hoses
87	K
86	Kidney dish
141	L
108	Labofill - cordless electronic pipette filling device labofill
64	Lining for glass tapers
115	Long spout dropping bottle screw caps
106	Long-stem funnels
135	L-shaped spreaders
137	"L" untapered connectors
18	M
18	Magnetic stirrer, non-electric
87	Male bed bottle with cap
71	Measuring jugs, long form
32	Measuring jugs, short form
70	Measuring scoops
66	Metal trays for histology
24	Micropipettes - pluripet ii
28	Microscope cover
71	Microscope slide boxes
97	Microscope slide storage and dust cover
50	Micro test tube rack preparation
43	Micro test tubes
54	Micro test tubes with caps for cobas-bio® and cobas mira®
107	Microtitre plates
107	Minifix
49	Mono-ball for vacuum with two valves
102	Mono-ball with valve and hoses
102	Multichannel variable volume micropipettes - multipet
36	N
37	Narrow neck bottles - grey
37	Narrow neck bottles with white tamper-evident cap
66	Non-return valve
20	O
20	Orange k32 hoses
24	Overturnable septa

H

INDICE ALFABETICO

ALPHABETICAL INDEX

INDEX ALPHABÉTIQUE

P	
Pasteur pipette	125
Pestles serrated	27
Pipette/burette rinsing set, automatic	74
Pipette filler with adapter	76
Pipette filler with plunger	77
Pipette jar	75
Pipette pumps	77
Pipette stand	75
Pipette stand, vertical	75
Pipette tray for drawer	76
Policemen stirring rods	29
Powder funnels	28
R	
Rectangular base	72
Rods for clot detachment and extraction	125
Round bottom flask	98
Round bottom flask holder	72
Round bowl	38-30
S	
Safety box	32-84
Safety box rack	32-84
Safety shields	23
Sample containers	97
Sample cup stoppers	104
Sample tube	98
Sample vials	26-60
Sampling bottle for urine	86-87
Schilling burettes	54
Scintillation vials	103
Screw cap container	108
Screw cap containers	26-60
Sedimentation imhoff cone	54
Semitransparent empty boxes for refills	119
Separating funnel holder	70
Seropet	140
Set of 7 adapters (guko)	24
Short spout dropping bottle screw caps	39
Slide dispenser	57
Slide mailer	57
Slide storing system	56
Snap-on lid containers	61
Snap-on lids containers	110
Spare "o" rings	65
Spare parts set and microchannel micropipettes stand	135
Spare vacuum retention valve	65
Specimen container	107
Spray bottle	40
Sputum collection container	108
Stackable tanks	31
Staining jar hellendhal type	58
Staining jar schifferdecker type	59
Staining jar with 2 lids	58
Staining rack	58-59
Standardized screw caps for bottles din 168	38-39
Standardized screw caps for bottles din 168	53
Sterile disposable pipette tips with filter	121
Sterilization indicator tape	33
Stirrer blades for balloon flasks	90
Stirring ball, magnetic	92
Stirring bars, crosshead	92
Stirring bars, cylindrical	91
Stirring bars, disc	93
Stirring bars, double ended	92
Stirring bars, magnetic	91
Stirring bars, magnetic with pivot ring	93
Stirring bars, oval	92
Stirring bars retrievers, magnetic	93
Stirring paddles	28
Stirring rods	28
Stopcocks	22
Stoppers for disposable test tubes	106
Stoppers with hoses connections	23
Storage box for 4 racks	59
Straight connectors tapered	17
Supple connectors	91
T	
Tanks with spigot	46
Tape	94
Tapered stoppers	94
Technicon®, beckman®, konelab® sample cup	104
Technotrate accessories	133
Technotrate digital burette	133
Test tube caps	81
Test tube racks	81
Test tube racks two/three tier	83
Test tube racks two-tier	83
Test tubes	94
Test tubes with screw caps	105
Test tubes with screw caps	109
Test tube trays	83
Timers	22
Timers	22
Tips, disposable for automatic micropipettes	111
Tissue embedding cassettes, uni type	127
Tissue embedding cassettes mega type	127
Tissue embedding sponge	127
Tissue grinders	27
Tongue depressor	86
Tray for microscope slides	58
Trays	30
Triangular stirring bars, magnetic	93
Tube-rack, microcentrifuge (1,5 ml)	82
Tubes, with rim	106
"T" untapered connectors	16
U	
Unequal straight connectors tapered	17
Universal safety pipette bulb	76
Universal test tube rack	82
Urine cups with screw cap	109
Uv range cuvettes	100
V	
Vacuum hoses k120	21
"Vadosa" vacuum pipetting device	77
Variable volume autoclavable	138
Variable volume autoclavable micropipettes - pluripet	134
Viscous liquid pumps	67
Visible range cuvettes	100
Volumetric flasks	97
Volumetric flasks with cap	52
Volumetric flasks with	
screw cap	52-53
W	
Wash bottles	40
Wash bottle screw caps	39
Watch glasses	27
Watch glasses	91
Water jet pump	65
Weighing bottles	26
Wide mouth wash bottles, graduated	41
Wide neck bottles	96
Wide neck bottles with black tamper-evident cap	37
Y	
"Y" untapered connectors	16

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HOSES CONNECTORS

PP

Autoclavable. Dimensions are shown as both nominal and exact as elasticity of hoses will influence the choice of size. Use hoses selector Art.1759 to determine "grip". This will then give equivalent crest dimension for the particular connector concerned. Available in packages of 10 or 100 pcs.

**TESTED ACCORDING TO REACH
REGULATION 1907/2006 ANNEX XVII.**

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"T" UNTAPERED CONNECTORS

PP

Excellent chemical and temperature resistance. With serrations to hold hoses tightly and prevent slipping. Autoclavable.

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"Y" UNTAPERED CONNECTORS

PP

Useful for any hoses system and vacuum line.

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UNEQUAL STRAIGHT
CONNECTORS - TAPERED

PP

Inlet and outlet have different diameters.

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DISCONNECTORS - TAPERED

PE

These quick disconnectors are useful for joining hoses of glass, rubber or plastic. Connectors fit tightly together. Replace expensive adapters.

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DRYING TUBES - TAPERED

PP

Ideal for drying gases. The two ends have push-on tapered connections to take rubber hoses of 8,9,10 mm diameters.

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FLOW INDICATOR, BALL

SAN

For liquid flow from 50 ml/min. in horizontal position, and 140 ml/min. in vertical position. Minimum air flow 3 litres/min. horizontal and 8 litres/min. vertical. Speed of rotation of ball proportional to flow. Not to be used with corrosive fluids or over +50°C.

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Art.	Dimensioni mm Dimensions mm Dimensions mm	Conness. O.D. mm Connect O.D. mm Connect. O.D. mm
538	102x14x57h	6,5 - 10

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FLOW INDICATOR, PADDLE WHEEL

SAN

For liquid flows down to 5 ml/sec. Wheel spins freely and speed of rotation is proportional to flow rate. Maximum recommended flow rate up to 95 ml/sec. Max 2 bar pressure. Not to be used with corrosive fluids or over +85°C. Can be used with gases.

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Art.	Dimensioni mm Dimensions mm Dimensions mm	Conness. O.D. mm Connect O.D. mm Connect. O.D. mm
539	90x15x40h	6 - 11

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MEASURING SCOOPS**PP**

PP scoops shaped conveniently for sampling. Flat bottom with beveled angles for easy cleaning.
SUITABLE FOR FOODSTUFF.

Art.	Capacità ml Capacity ml Capacité ml	Dimensioni mm Dimensions mm Dimensions mm
1057	10	32x103x16
1058	25	43,8x137,3x21
1059	50	55,4x163,7x26
1060	100	67,2x201,7x34
1062	250	94x261,5x43,7
1063	500	112x315x55
1064	1000	141x386,6x69,4

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INDUSTRIAL SCOOPS**HDPE**

This scoop, made of HDPE, is particularly designed for heavy industrial use. Features a round shaped design to facilitate complete transfer of sampling material. **SUITABLE FOR FOODSTUFF.**

Art.	Capacità ml Capacity ml Capacité ml	Lunghezza cm Length cm Longueur cm
1110	250	25
1111	500	30
1066	1000	35
1112	1500	40

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ISOFLEX LABORATORY HOSES**PVC NON-TOXIC, AUTOCLAVABLE**

Autoclavable PVC hoses, transparent non-toxic; non ageing and non oxidizing. Supplied in 20 meter rolls packed in PE bags. The hoses are marked every meter for ease of use.

Technical data

Colour: clear
Odour: none
Taste: none
Toxicity: non-toxic
Ageing: none
Oxidation: none

Chemical resistance

Strong acids: good
Weak acids: excellent
Strong alkalis: good
Weak alkalis: excellent
Organic solvents: fair

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HOSES K30

PVC

Flexible and transparent hoses manufactured in atoxic material. Suitable as components for medical devices, industrial and hospital products. DEHP free. Temperature: -15°C/+60°C. Conform to PhEur 3.1.1.1 / USP XXXII 2009. **SH. 65+/- 5**

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HOSES K70

SYLICON

Flexible and transparent, suitable as components for medical devices. Conform to FDA 21CFR§177,2600, PhEur 3.1.9, ISO 10993-5. Temperature: -60°C/+200°C. **SH A 60 +/- 5**.

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HOSES

PTFE

Excellent resistance to acids, bases and solvents. Temperature up to +260°C. Autoclavable. **SH D 55/65**.

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ORANGE K32 HOSES

NATURAL RUBBER

Utilful for gases, liquids and medium concentration acids. Temperature: +70°C/-35°C. **SH A 60 +/- 5**.

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VACUUM HOSES K120**PARA RUBBER**

Hoses with cloth impression surface. Suitable for air, water, medium concentration acids and alkaline solutions, abrasive products, oxygenated solvents etc. Temperature: +70°C/-40°C. SH 45 +/- 5.

Art.	Dim. mm	PKG. m
3885	5x15	6
3886	6x18	6
3887	8x16	6
3888	12x21	6

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HOSES**NEOPRENE**

For normal laboratory use. Resistant to acids and solvents, benzole and similar excluded. Temperature: +85°C/-35°C. SH A 65 +/- 5.

Art.	Dim. mm	PKG. m
3880	4x6	20
3881	6x9	20
3882	8x12	10
3883	10x14	10

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HOSES**LATEX**

Hoses with smooth and transparent surface. Density 0.92. Autoclavable. SH A 40 +/- 5.

Art.	Dim. mm	PKG. m
3870	4x7	10
3871	6x9	10
3872	8x12	10
3873	12x17	10

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Permette la selezione dei tubi in plastica, vetro o gomma secondo il diametro interno ed esterno.
Diametri int. disponibili: mm 1 / 2,5 / 3 / 4 / 4,5 / 5 / 6 / 7 / 8 / 9,5 / 10 / 11 / 12 / 13 / 14 / 15 / 16 / 17 / 18 / 19 / 20 / 21 / 22 / 25.
Diametri est. disponibili: mm 1,5 / 2,5 / 3 / 4 / 5,5 / 6 / 7 / 8 / 9 / 10 / 11 / 12 / 13 / 14 / 15 / 16 / 17 / 18 / 21.

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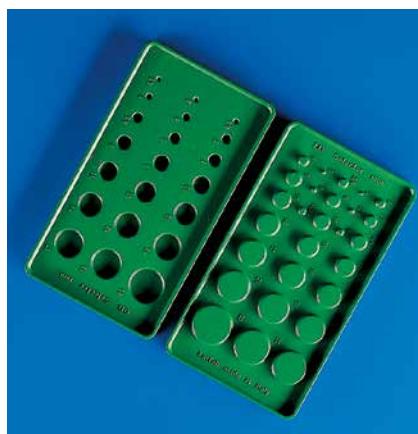
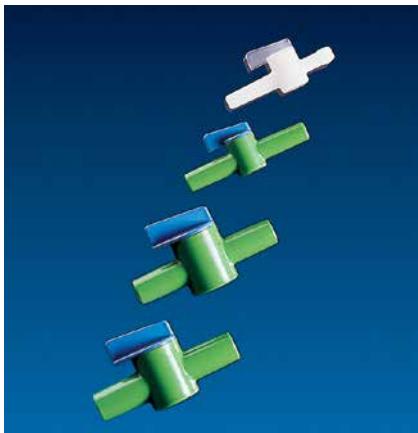
STOPCOCKS

PP AND HDPE

Ideal for liquids and gases under low pressure (operating pressure 0,5 bar) or low vacuum. HDPE body interfaces with PP tap giving a smooth turning action requiring no lubricants. All are straight bored.

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TIMERS

ABS CASE

The Clip model is portable and pocket size, with mechanical movement (accuracy 3%) and timing up to 60 minutes. The Onehand has the same performance as the Clip but with the advantage that it can be rewound using one hand only.

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ELECTRONIC TIMERS

ABS CASE

It is battery operated and has a 4 digits display up to 19 hours 59 minutes. An alarm sounds at the end of the set time. Supplied with support, clip and magnet.

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SAFETY SHIELDS

PMMA

The safety shields are available in two sizes with different thicknesses and are made of high quality, clear acrylic; are free standing, portable and stable, and are all fitted with non-skid feet. Art. 571 is designed to provide laboratory and healthcare personnel with facial protection from hazardous substances, aerosols and splashing. Art. 572 is designed to provide additional protection from radiations while using beta emitting isotopes, including P32.

Art.	Dim. mm	Spessore mm / Thickness mm / Epaisseur mm
571	242x365h	5
572	305x480h	9

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STOPPERS WITH HOSES CONNECTIONS

PP

Standard taper 1:10 ensuring perfect sealing action. Hose-adapters Ø mm 8 - Length 20 mm.

Art.	Cono NS NS Cone Cône NS	Ø inf. mm lower Ø mm Ø inf. mm
1816	29 / 32	26,2

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CONICAL STOPPERS

RED RUBBER

General-purpose laboratory red rubber stoppers. Autoclavable and pierceable

Art.	Ø sup. mm upper Ø mm Ø sup mm	Ø inf. mm lower Ø mm Ø inf. mm	h mm	Art.	sup. upper Ø mm	inf. lower Ø mm	h mm
1129	9	6	18	1139	30	21	29
1131	12	9	18	1140	32	25	34
1132	13	10	18	1142	37	28	37
1133	15	11	18	1143	42	32	41
1134	16	12	20	1144	45	34	45
1135	18	13	22	1145	48	37	50
1136	20	14	24	1146	54	41	53
1137	23	16	26	1147	60	43	60
1138	26	19	28	1148	65	52	70

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CONICAL STOPPERS

SYLICON

Autoclavable.
SUITEABLE FOR FOODSTUFF.

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OVERTURNABLE SEPTA

SYLICON

Pierceable stoppers to withdraw solutions and suspensions from closed containers, using syringes. Suitable for use in food and beverage industry.
H: height of overturnable section
h: height of lower section
SUITABLE FOR FOODSTUFF.

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GAS SAMPLING BALLOONS

LATEX

Latex balloons with volume range from 1,000 cc to 8,000 cc

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**MONO-BALL WITH VALVE AND
HOSES****RED RUBBER**

Suitable for Pellet burette charging or for vaporizers.

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DOUBLE BALL WITH NET**RED RUBBER**

Suitable for automatic burettes and flash chromatography columns.

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DROPPING BULBS WITH RIM**LATEX**

Suitable for Ranzier dropping tubes and Pasteur pipettes.

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HOT GRIP

SYLICON RUBBER

It provides protection against hot containers up to +260°C and cold containers down to -57°C. Internal relief surface.

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SAMPLE VIALS

LDPE

General purpose containers with push on hinged lids. Flat bottomed. Lightweight and unbreakable. Can be used as weighing containers.

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SCREW CAP CONTAINERS

LDPE

Ideal for storing mercury etc. Leakproof. Small quantities of liquid or solid samples can be taken out by pipette/spatulas without the hindrance of internal shoulders. Sturdy, heavy walled with excellent chemical resistance.

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WEIGHING BOTTLES

PP

Reduce the tare weight correction with these light weight, autoclavable, reusable translucent bottles. Airtight push on lids included. Highly resistant to acid attack. No internal shoulders to allow use of spatulas. Tops can be squeezed to assist pouring. Remove lid before autoclaving.

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WATCH GLASSES

PP

Very stable having a moulded base. Can withstand temperatures up to 120°C.

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TISSUE GRINDERS

BOROSILICATE GLASS

Tissue grinder is composed by two components to be ordered separately: Borosilicate glass mortar (items 6102 to 6150) and PTFE pestle (items 6302 to 6350). Ideal for small amounts of tissue. PTFE pestle has 0.15 - 0.25 mm clearance with borosilicate mortar. Volumes stated are working volumes. Avoid excessive generated heat as could cause binding.

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PESTLES SERRATED

HEAD MADE OF PTFE, SHAFT MADE
OF STAINLESS STEEL

Pestle should be used in combination with borosilicate mortar for sample crushing. Ideal for tough materials such as tumors. Pestle shaft diameter is 6,5 mm.

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MAGNETIC STIRRER,
NON-ELECTRIC

PP - ABS

This unique stirrer is operated either by water pressure between 1 and 2.5 Kg/cm2 (15 to 40 psi) or by air pressure. No electric supply is needed thus eliminating heat, shock and explosion hazard. Corrosion proof allplastic construction. Can be used on the bench or clamped to a stand with 12 mm (1/2") diameter (our part. 267). Supplied complete of tubes.

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POWDER FUNNELS

PP

60° angle, autoclavable funnels. Large bore avoids bridging of powders during transfer. Ribbed outside to avoid pressure lock. Can be used with liquids. Excellent chemical resistance.

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GENERAL PURPOSE FUNNELS

HDPE

Tough, robust, general purpose funnels suitable for a wide range of liquids. Upstand on rim to reduce possible overflow. Ribbed on outside to eliminate back pressure. Hanging eye for storage.

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STIRRING PADDLES

PP

The blades are made of chemically resistant PP and are fixed to a steel rod covered with PP with a diameter of 8 mm x 350 mm long.

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STIRRING RODS

PVC

Rigid, light and unbreakable. The ideal replacement for glass rods.

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POLICEMEN STIRRING RODS

PP

White, one-piece rod with flattened ends forming two small spatulas.

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ICE BUCKET

EXPANDED PUR

Equipped with lid and two side-handles. Thickness of about 15 mm. Red colour. Excellent isolating feature. Can be used with dry or minced ice and cooling mixtures. Liquid nitrogen resistant for a few minutes.

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INPUT TRAYS

PVC

Very useful for knick knacks - rulers, erasers, pencils etc, in desk drawers. Central compartment of five compartment tray can be used for pipettes. Sturdy. Can be used for carrying items such as stir bars, connectors etc.

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INPUT TRAYS

PP/PVC

Art. 548 is a tough 12 liters autoclavable tray for a variety of uses. Can hold 18x500 ml or 12x1 lt bottles.

Art. 952 is a strong PVC tray for holding sample bottles with base diameter not greater than 22 mm. Ideal for 7 ml (1/4oz) Bijou bottles or specimen tubes. Each of nine rows can hold 12 of such bottles making a total capacity of 126.

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DEEP TRAYS

PVC

With the ribbed bases, these are ideal for photographic purposes. Stackable.

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ROUND BOWL

PP

White bowl with rim, autoclavable at 121°C for 20 minutes. Sturdy structure. Two grips have been designed in the lower part of the rim to help carriage.

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TRAYS

HIGH IMPACT PS

Suitable for foodstuffs and general laboratory use.

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DEEP TRAY, LOW FORM, STACKABLE

HDPE

Stackable deep trays made of HDPE with handles.

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**DEEP TRAY, LOW FORM WITH
HANDLES AND DRAINING HOLES****HDPE**

Rectangular tank with draining holes on all sides. Stackable with recessed rim to allow ventilation.

Art.	Dim. mm	h mm	Cap. It
547	350x540	115	16

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STACKABLE TANKS**HDPE**

Can be stacked. High resistance to organic and inorganic acids. Robust with lifting lugs and strengthening ribs to ensure long life.

Art.	Dim. ext. mm	h mm	Cap. It
602	315x450	200	20
603	370x625	255	46
604	440x730	310	72
605	500x600	445	100

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GREY TANKS**HDPE**

Useful general purpose tanks that can be stacked for storage.

Art.	Cap. It	Dim. mm	Colore / Colour / Couleur
3371	27	340x455x250	Grigio / Grey / Gris
3372	40	350x560x304	Grigio / Grey / Gris
3373	75	460x760x310	Grigio / Grey / Gris

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BUCKET WITH SPOUT**LDPE**

Graduated bucket by liter with pouring spout.

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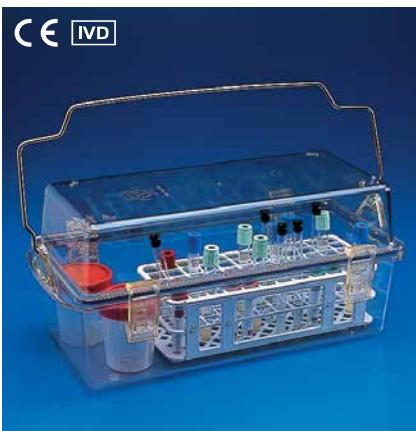
SAFETY BOX

PC

Made of polycarbonate and fully autoclavable. Designed for transporting tubes, containers, vials, plastic bags and ready-to-use syringes whenever an airtight seal is required. Designed to provide a high degree of safety when transporting antineoplastic drugs from the Preparation Center to the site where they will be administered. Completely clear to enable inspection for possible container leakage before opening. The cover has a hermetic silicon safety seal to ensure containment of fluids or anti-tumor drugs and protects against leakage and/or accidental spill from bottles or seepage from plastic bags. Closes using 4 safety locks preventing accidental opening if the box is dropped. Inside the box, beveled edges and the special design of the bottom guarantee safe, simple and efficient cleaning. An indelible biological hazard symbol is moulded on the outside of the cover. Includes a stainless steel AISI 304 handle for ease of transport and an instruction sheet on use and maintenance.

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SAFETY BOX RACK

STAINLESS STEEL

Designed for 10 urine containers up to 200 ml size and 4 stool containers 30 ml size.

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FRIGO BOX, SQUARE

HIGH IMPACT PS

Low temperature containers made of high impact thermoplastic material, thermally isolated by means of a polyurethane isolating jacket; non-toxic and therefore 100% foodsafe. A 1 Kg cooling block, art. 3707, is available for art. 3705, 3706. Inside art.3706 1pc. art. 569.

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STERILIZATION INDICATOR TAPE

ADHESIVE PAPER TAPE

Art. 1052 is made by impregnated semi-creped paper and adhesive natural rubber. To be used with individual units (i.e. packs, containers) to confirm sterilization treatment and to distinguish between processed and unprocessed units. After sterilization, indicator strips change from clear colour to brown. (minimum conditions for change are: steam 10 min 121°C or 2 min 134°C). Lead free. If correctly stocked, it can be used till 3 years from production date.
NOT INTENDED FOR USE AS VALIDATION SYSTEM OF STERILIZATION PROCESS.

Art. 1051 is an adhesive tape with good resistance to heat and damp; it's particularly recommended for indicating dry sterilization (POUPINEL method). Made by semi-creped, impregnated paper and adhesive Silicone it is easy to remove after treatment. After sterilization the "V" indicators strips changes colour from light green to brown (minimum conditions for change are: dry heat 30 minutes at 155°C). Lead free. If correctly stocked, it can be used till 2 years from production date.
NOT INTENDED FOR USE AS VALIDATION SYSTEM OF STERILIZATION PROCESS.

Both Tapes are in conformity with ISO 11140/2005

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GRADUATED WIDE NECK BOTTLES

PE (BODY) PP (CAP)

Will withstand considerable pressure before rupture. Secured by tagging points on bottle and cap. Confidential notes can be placed in the space between insert sealing plug and cap.

SUITABLE FOR FOODSTUFF according to:

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GRADUATED SQUARE BOTTLES WIDE NECK

PE (BODY) PP (CAP)

Space saving bottles, ideal for long term storage of samples. Approximate graduations allow reproducible filling without the use of cylinders/beakers. Easy grip caps can be wired down for security. **SUITABLE FOR FOODSTUFF** according to:

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GRADUATED NARROW NECK BOTTLES

PE (BODY) PP (CAP)

Double sealing with insert cap to assure tightness. Sloping shoulders reduce tendency for residue deposits. No leaking. Tagging points on shoulder and cap ensures tamper proof closure. Easy grip cap.

SUITABLE FOR FOODSTUFF according to:

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NARROW NECK BOTTLES - GREY**PE**

Completely opaque. Excellent for all photosensitive material whether liquid or solid. Sloping shoulders reduce tendency for residue deposits. No leaking. Insert plug for maximum security. Easy grip cap.

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**NARROW NECK BOTTLES WITH
WHITE TAMPER-EVIDENT CAP****HDPE**

The tamper-evident cap incorporates a safety ring which is secured when the cap is screwed on to the bottle. Opening the bottle breaks the safety ring, showing signs of interference.

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**WIDE NECK BOTTLES WITH
BLACK TAMPER-EVIDENT CAP****HDPE**

The tamper-evident cap incorporates a safety ring which is secured when the cap is screwed on to the bottle. Opening the bottle breaks the safety ring, showing signs of interference.

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GRADUATED WIDE NECK BOTTLES

PP

Autoclavable, graduated bottle meeting Food and Drug regulations with pictograms. Very sturdy, ruptureproof bottle. Long neck thread ensures leakproof closure. Wide neck for easy filling and emptying of liquid or powder samples. Conforms to DIN 13316 and 168.

SUITABLE FOR FOODSTUFF.
according to:

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GRADUATED NARROW NECK BOTTLES

PP

Autoclavable, graduated, heavy walled bottle. Ideal for transporting toxic waste waters, boiler feed water etc., back to laboratory for analysis as samples can be secured using shoulder and cap tag to avoid sample contamination in transit. Conforms to DIN 13316 and 168. Remove cap before autoclaving.

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STANDARDIZED SCREW CAPS FOR BOTTLES DIN 168

PP

Standardized screw caps for PP bottles DIN 168. Cap tag to avoid sample contamination in transit.

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GRADUATED NARROW NECK BOTTLES - WITHOUT CAPS

PE

An ideal bottle for water and oils - there are no shoulders to restrict the flow. They are graduated and so can be used as rough measures. Can be secured using shoulder and cap tag to avoid sample contamination in transit. Conforms to DIN 13316 and 168.

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**SHORT SPOUT DROPPING BOTTLE
SCREW CAPS****PE**

They fit narrow neck standardized bottles part. no. 15801 - 15821 - 15841 - 15861 - 15881 - 15901 and wide neck standardized bottles part. no. 1608 - 1610. Supplied with PE cap for tip closure.

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GL25

GL32

Lungh. beccuccio/Spout length/Long. bec mm

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DISPENSING BOTTLE

HDPE-PP

Suitable for liquids single and/or continuous dispensing.

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WASH BOTTLES

PE

Ideal for distilled water, the stream of liquid is regulated by hand pressure. Dispensing tip gives ultra-fine stream or can be removed to increase flow.

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WASH BOTTLES

PE

The stream of liquid is obtained by blowing through the straight tube. The direction of the stream of liquid can be controlled by hand.

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**DROPPING BOTTLES -
RANVIER TYPE**
PE

Complete with a PE tube with rubber teat.
SUITABLE FOR FOODSTUFF.

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ADATTE AL CONTATTO ALIMENTARE.

INTEGRAL WASH BOTTLES
PE

Easy squeeze wash bottles with integral moulded tube. Dispensing tip, which is provided with PE closure cap, can be cut back to increase flow. Cap can be use also with bottles series 15801 - 1631. Special shape for easy grip and stability.

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ADATTE AL CONTATTO ALIMENTARE.

**WIDE MOUTH WASH BOTTLES,
GRADUATED**
PE

Wide mouth for easy, safe filling. Colour coded caps for easy identification. Graduated. Ingenious cap designed to avoid involuntary dripping due to pressure build up in bottle - where toxic or radioactive material is being used. Available in the colours: 00 neutral, 04 blue, 06 yellow, 10 red. For example: 1637-10 is a red capped 250 ml wash bottle.

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ADHESIVE LABELS

Transparent, punched labels for use with wash bottles and plastic or glass reagent bottles, as recommended by the NFPA (National Fire Protection Agency). Complete with risk assessment numbers, recommendations for use and special notices shown in reagent catalogues.

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CYLINDRICAL JARS WITH RIBBED CAP

PE - JAR AND INSERT
PP - CAP

The cylindrical jars are supplied with insert and cap. Thanks to the large neck they are suitable for powders, solid specimen, creamy or semisolid substances, crystals and hygroscopic substances.

SUITABLE FOR FOODSTUFF
according to:
 • EC Regulation 10/2011

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ADATTE AL CONTATTO ALIMENTARE.

**ADJUSTABLE VOLUME
DISPENSER BOTTLES****PE (BOTTLE)
PMP (CONTAINER)**

The graduated clear PMP measuring cup slides up and down the tube and the variation in height changes the volume to be dispensed. The dispensed volume remains constant at any preset position of the cup. The cup is filled by squeezing the bottle and forcing the liquid up the tube. When the bottle pressure is released, the excess will be drawn back into the bottle and only the desired volume will remain.

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GRADUATED BOTTLE**HDPE**

Specific design with rectangular base thought for easy storage of the bottles on the shelves and for enhanced stability when in use; the ergonomic design of the handgrip allows easy pouring. Screw capped and graduated up to 2.000 ml.

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BOTTLE-CARRIERS**PP**

Designed to carry 6 and 4 bottles, with maximum diameter up to 90 mm and 110 mm respectively. Both of them, made of PP, feature an ergonomic design of the handle and can be easily stacked.

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FIRST EYD EMERGENCY EYE WASH STATION

HIGH IMPACT PS

Emergency eye wash station moulded in strong high-impact PS. The directions are clearly printed on the support rack and the bottle can be filled with distilled water or other medicated liquid for ophthalmic use. Eye wash bottle art. 383, to be ordered separately, is made of PE provided with an eye cup with snap on cap, carefully designed to allow a quick and immediate flushing to remove foreign matters and prevent serious eye damage. The head has an air-vent valve to assure air refill of free space in the bottle thus preventing draw back of contaminated wash into the bottle. A perforated spray plate in the cup centre breaks up the stream to give the eye a gentle wash. A drain tube releases waste.

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INTEGRAL EYE WASH BOTTLE M45

PE

Can be filled with distilled water or any other medicated liquid intended for ophthalmic use. The bottle has an eyecup with a specially designed snap-on cap for fast and convenient eye washing to remove foreign objects and prevent serious damage to vision. An automatic valve is incorporated in the head to admit air and prevent the contaminated eyewash from flowing back into the bottle. The eyecup has a perforated plate to break up the stream and provide a uniform spray to the eye. Also includes a tube to drain the contaminated eyewash.
SUITABLE FOR FOODSTUFF.

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**HEAVY-WALLED CARBOY
BOTTLES, WIDE NECK****HDPE**

Very resistant bottles, ideal for liquid or powders. Wide mouth for easy filling and emptying. Insert plug added. Caps can be wired down for safety. Tommy bar lugs on 10 litre capacity to assist release. Carrying handles.

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**HEAVY-WALLED CARBOY
BOTTLES, NARROW NECK****HDPE**

Very resistant heavy-walled carboy bottles. O-ring seal cap make them leakproof. The cap can be wired down for added safety. Complete with carrying handles.

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**ASPIRATOR BOTTLES, HEAVY
DUTY****HDPE**

Designed for life time service. Can be used in sub zero temperatures. Very sturdy having thick walls whilst retaining translucency. The 3/4" BSPT spigot (supplied) mates with a 12 mm deep thread in the aspirator for total leakproof performance. Recommended that a few drops of vaseline be put on the thread before assembly.

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TANKS WITH SPIGOT

HDPE

These two products, with size 10 liters and 20 liters, are specifically designed to rationalize laboratory and small production line organization, giving the opportunity to optimize water, aqueous solutions and solvents availability in a reduced space. They are made of High Density PE food grade, with a comfortable handle for displacement. The spigot is placed in a position allowing the tank to stand, as to facilitate the use of these tanks in packaging and transporting liquids.

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INDUSTRIAL TANKS

HDPE

These four tanks feature a standardised industrial packaging design, with sizes 2, 5, 10 and 20 litres, and even if specifically addressed to a use in industrial packaging, they could result useful in everyday laboratory practice. They are made of High Density PE food grade, with a comfortable handle, and feature a secure seal-cap following DIN recommendations. They are conform to UN regulation on transportation of dangerous and toxic liquids and are suitable for road (ADR-RID), sea (IMO-IMDG) and air (ATA-ICAO) transport.

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GRADUATI E TARATI

GRADUATED AND VOLUMETRIC PLASTICWARE

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GRADUATED TALL FORM MEASURING CYLINDERS - CLASS B

PP

Conforms to ISO 6706 - 1981 (E). Special PP giving very high translucency. Autoclavable to 121°C for 20 min., continuous work at 100°C. Permanent graduations are moulded in for reproducibility. No wetting. Chemically cleaner than glass. Chemically non-absorbant. Excellent chemical resistance. Pentagon base for added stability.

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GRADUATED TALL FORM MEASURING CYLINDERS - CLASS B

PMP (TPX°)

Conforms to ISO 6706 - 1981 (E). Crystal clear. Autoclavable. Can be used for liquids up 170°C. No wetting. No meniscus. Permanent graduations are moulded in for reproducibility. Chemically cleaner than glass. Chemically non-absorbant. Excellent chemical resistance. Pentagon base for added stability.

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BLUE GRADUATED MEASURING CYLINDERS TALL FORM - CLASS B

PP

Conforms to ISO 6706 - 1981 (E). Special PP giving very high translucency. Autoclavable to 121°C for 20 min., continuous work at 100°C. Permanent graduations.

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BLUE GRADUATED MEASURING CYLINDERS TALL FORM - CLASS B
PMP (TPX®)

Conforms to ISO 6706 - 1981 (E). Crystal clear. Autoclavable. Can be used for liquids up 170°C. No wetting. No meniscus. Permanent graduations. **SUITABLE FOR FOODSTUFF.**

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GRADUATED MEASURING CYLINDERS, SHORT FORM
PP

Low form, circular base. Excellent chemical resistance. Autoclavable. Permanent graduations moulded into cylinders. Can withstand temperatures up to 120°C.

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GRADUATED MEASURING CYLINDERS, SHORT FORM
PMP (TPX®)

Low form, circular base. Excellent chemical resistance. Autoclavable. Permanent graduations moulded into cylinders. Can withstand temperatures up to 170°C.

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GRADUATED BEAKERS,
LOW FORM - CLASS B

PP

Conforms to ISO 7056 - 1981 (E). New PP giving very high translucency. Autoclavable to 121°C for 20 mins. continuous work at 100°C. Permanent graduations are moulded in for reproducibility. Excellent chemical resistance.

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GRADUATED BEAKERS,
LOW FORM - CLASS B

PMP (TPX®)

Conforms to ISO 7056 - 1981 (E). Crystal clear, autoclavable. Will withstand 170°C for short periods. Permanent graduations are moulded in for reproducibility. Excellent chemical resistance.

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BLUE GRADUATED BEAKERS,
LOW FORM - CLASS B

PP

Conforms to ISO 7056 - 1981 (E). New PP giving very high translucency. Autoclavable to 121°C for 20 min., continuous work at 100°C. Permanent graduations. Excellent chemical resistance.

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BLUE GRADUATED BEAKERS,
LOW FORM - CLASS B

PMP (TPX®)

Conforms to ISO 7056 - 1981 (E). Crystal clear. Autoclavable. Will withstand 170°C for short periods. Permanent graduations. Excellent chemical resistance.

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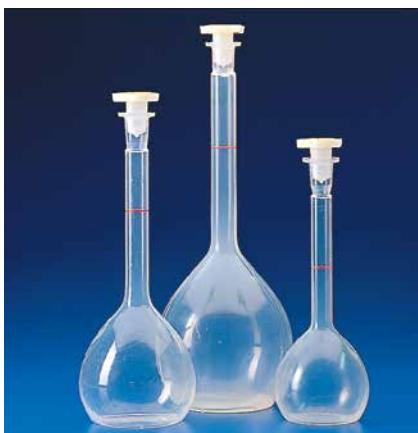
VOLUMETRIC FLASKS WITH CAP

PP

These flasks create no meniscus so final readings are easy. Calibrated gravimetrically at 20°C. Translucency of PP enhanced when in contact with liquid. Thick walls eliminate distortion. Thin neck make final reading precise. Autoclavable. Chemically non-absorbant. No wetting. Stoppers included.

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VOLUMETRIC FLASKS WITH CAP

PMP (TPX®)

These crystal clear volumetric flasks create no meniscus so final readings are easy. Calibrated gravimetrically at 20°C. Thick walls eliminate distortion. Thin necks make final measurement precise. Autoclavable to 170°C. Chemically cleaner than glass and chemically non-absorbant. No wetting. Stopper included.

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CONICAL STOPPERS

PE

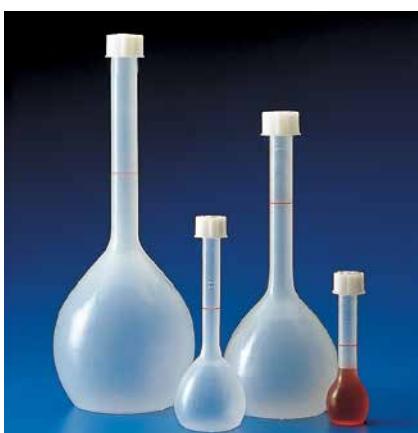
Stoppers manufactured to DIN Standards 12242 ensuring sealing action on non-standard surfaces.

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VOLUMETRIC FLASKS WITH SCREW CAP

PP

These flasks create nomeniscus so final readings are easy. Calibrated gravimetrically at 20°C. Translucency of PP enhanced when in contact with liquid. Thick walls eliminate distortion. Thin neck make final reading precise. Autoclavable. Chemically non absorbant. No wetting. Screw caps included.

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GRADUATED, SEPARATING FUNNEL**PMP (TPX®)**

Crystal clear funnel with permanent graduations moulded into body. Needle control for precise control of flow (important for end point). No wetting. Top unscrews and valve can be detached for thorough cleaning. Autoclavable. Excellent chemical resistance. Complete with stopper.

Art.	Cap. ml	Grad. ml	to level ml	NS DIN
978	500	2,0 5,0 25,0	20 50 500	29 / 32

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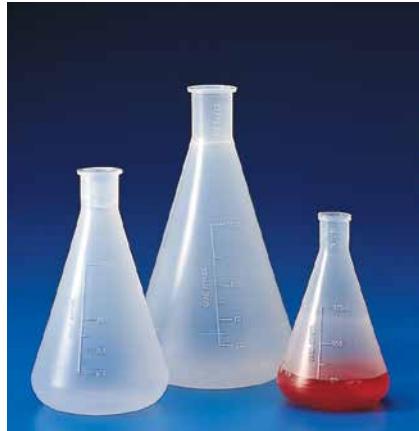
SEDIMENTATION IMHOFF CONE

SAN

Crystal clear Imhoff cone. Permanently moulded graduations. Easy draining via PP screw plug in the base. Back flush cleaning easily effected.

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SCHILLING BURETTES

PE / PMMA

Automatic plastic Schilling burettes (according to din 12.700 standards, Class B). With push button for drop-by-drop delivery. PE reservoir with stable PE-LD base.

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GRADUATED BURETTES

PVC / PMP / FEP

High precision plastic burettes (according to DIN 12.700 standards - Class B), particularly useful for routine work to eliminate the danger of glass breakages. The combination of different plastic materials (PMP, FEP, PVC) provides good resistance to chemicals. The stopcock assembly is moulded in one piece, with FEP key; requiring no lubricant and providing a leak-proof seal.

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MICROSCOPE SLIDE STORAGE
AND DUST COVER

ABS / PS

Ideal for storage of large quantities of 76x26 mm microscope slides. Each unit will accommodate 100 slides. Each cavity is numbered. Batch data can be written on either the inside or outside of the end plates. It is strongly recommended that the covers (Art. 920) are used to avoid dust and other foreign bodies. These can then be stored in the cabinets (Art. 930).

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MICROSCOPE SLIDE BOXES

HIGH IMPACT PS

The slots are well spaced for easy handling of 76x26 mm slides. Each slot is clearly numbered and can be cross referenced with corresponding index card. Lids are totally removable. The 50 and 100 place boxes have locating lugs so boxes can be stacked. Virtually unbreakable.

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SLIDE MAILER

PE

Strong protection for 76x26 mm specimen slides during transportation. Each slide has its own compartment with support bars at 55 mm centres to protect subject matter from being scuffed. Particularly useful for cytologic and geological specimens.

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SLIDE MAILER

PP

Ideal for storing and shipping 5 to 10 microscope slides. Unbreakable, chemically inert, the container will also hold up to 12 ml as a specimen vial; 20 mm of slide projects above jar, allowing easy removal. The thread of cap is very deep allowing leakproof seal for shipment. Supplied with small loop on cap for tagging. Cap should be taped for shipment.

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SLIDE DISPENSER

ABS / PS

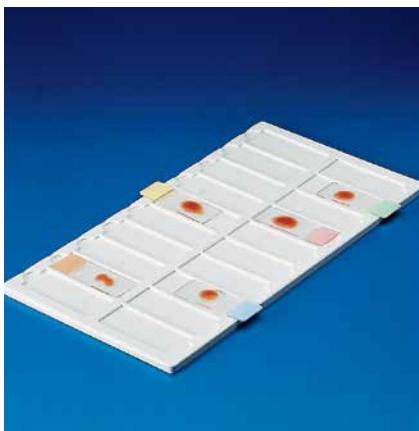
For finger print free dispensing of 76x26 mm microscope slides. Up to 50 slides can be put in the magazine. A cover protects the slides from dust. Each slide is dispensed by rotating the knurled knob on either side of the moulded base. Rubber rollers at the sides of the slides ease out the slide which can then be handled by the edges.

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TRAY FOR MICROSCOPE SLIDES

PVC

Moulded in white PVC for easy identification of different dyes on the slides. Ideal work station for 76x26 mm slides. Designed for easy removal of slides.

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STAINING JAR WITH 2 LIDS

PMP (TPX®)

Moulded in PMP for use with Art. 354. Supplied with 2 lids - one lid for use when dish is not in use, providing evaporation-free storage of staining liquid - the other lid with special opening to allow handle of staining rack Art 354 to be inserted; Xylene can cause etching after 2 days at room temperature.

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STAINING RACK

PP

Staining rack made of PP with horizontal tray, holds 20 slides 76x26 mm (3"x1"). For use with Art. 353.

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STAINING JAR HELLENDHAL
TYPE

PMP (TPX®)

PMP staining jar holds 76x26 mm (3"x1") slides vertically. Eight grooves. Supplied with PMP cover. Can accommodate 16 slides back to back. Xylene can cause etching after 2 days at room temperature.

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**STAINING JAR
SCHIFFERDECKER TYPE****PMP (TPX®)**

Staining jar for 76x26 mm (3"x1") slides. Ten grooves. Supplied with PMP cover. Can accommodate 20 slides back to back. Xylene can cause etching after 2 days at room temperature.

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STAINING RACK**POM**

Staining rack for simultaneous staining of 25 slides; to fit trough Art. 1101.

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25 SLIDES STAINING TROUGH**POM**

Staining trough for 25 slides; accomodating one rack, Art 1100. It can also be used as a mailing box.

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STORAGE BOX FOR 4 RACKS**PS**

Storage box only; can take up to 4 staining racks. Art. 1100, for a total quantity of 100 slides.

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SNAP-ON LID CONTAINERS

PS

Cylindrical containers in PS material, wide mouth with snap on lid.

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COLORIMETRIC TRAY

PE

With three round depressions. Particularly useful for colorimetric tests.

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MICROSCOPE COVER

PMMA

Clear acrylic resin microscope cover with a square base; protects against dust and dirt.

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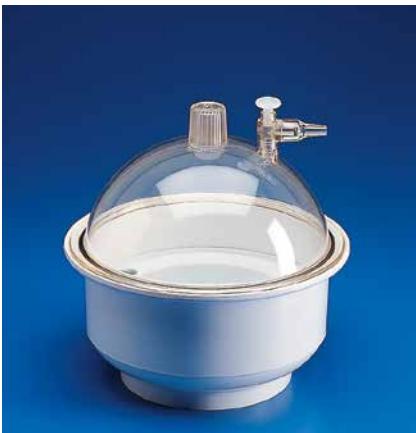
FILTRATION AND VACUUM DESICCATORS

PP / PC

Implosion proof desiccators. Will take vacuum of -740 mm Hg with loss (absolute zero: -760 mm/Hg) not exceeding 20 mm Hg in 24 hours. The transparent polycarbonate cover seals onto a neoprene 'O' ring set into the flange of the PP base. A little grease spread on the 'O' ring will assist in establishing initial vacuum. Within the desiccator is a removable pan for holding a drying agent. Vacuum release tap/non return valve assembly is included (Art. 229); specially grooved stopper in this assembly allows air, when required, to be re-admitted very slowly to avoid disturbing the contents.

Usage Instructions:

- remove the cover in PC - insert the samples in the pan
- open the vacuum retention valve (PE needle) connect the tube of the pump to the desiccator spigot in order to create vacuum close the vacuum retention valve

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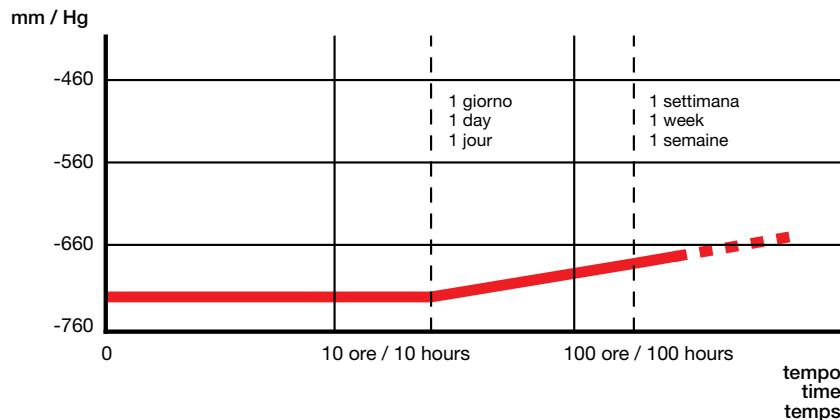
DESICCATOR PLATES

PP

These plates made of PP are extremely useful as a support for crucibles, Petri dishes etc. Use at room temperature. Not suitable for use with hot crucible.

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SPARE VACUUM RETENTION VALVE

PC WITH PE NEEDLE

Vacuum retention valve for desiccators
codes 550-230-554

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WATER JET PUMP

PP

This pump will work on pressures up to 10 Kg/cm². Built in non return valve to eliminate back flow. Useful for sucking to waste any spilt chemicals as well as for filter work. Easily dismantled for cleaning. Autoclavable.

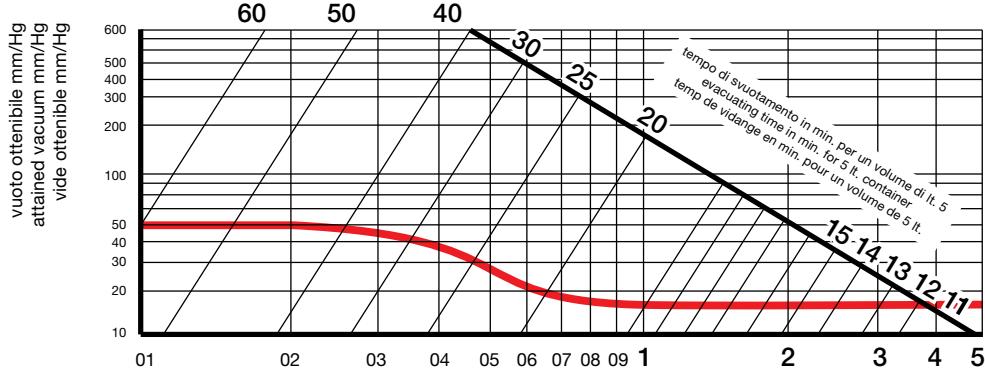
The graph below details the performance of the pump under different conditions. Example: Operating pressure of water: 1 Kg/cm². Attained Vacuum: 15 mm/Hg. Will evacuate a 5 litre container in about 20 minutes.

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NON-RETURN VALVE

PP INNER DISK: DUTRAL

Prevents water back-flow when pressure falls. Useful for any tubing system and vacuum line. Maximum safe working pressure: 2 BAR (equal approx. 28.4 P.S.I.). Minimum safe working pressure 0,07 BAR at room temperature (20°C).

Tested according to Reach Regulation 1907/2006 Annex XVII

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FUNNELS, BUCHNER

PP

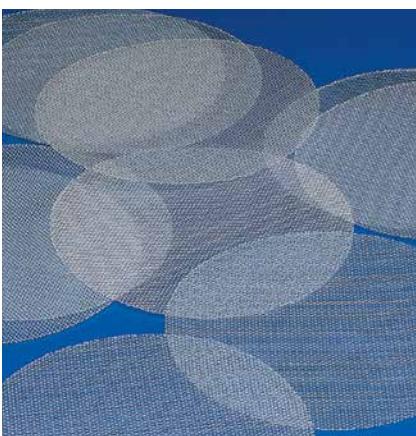
Robust, lightweight. Unbreakable and resistant to corrosion and heat. Two piece construction. Top and bottom can be separated for easy and total cleaning. Autoclavable. Excellent chemical resistance. SUITABLE FOR FOODSTUFF.

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DISKS FOR BUCHNER FUNNELS

PE

Useful interface between filter paper and perforated Buchner plate. Avoids clogging of filter holes. Texture 300 micron (μm).

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HAND OPERATED VACUUM PUMP

HIGH IMPACT PS

Light, portable hand operated vacuum pump. Attains a vacuum of 625 mm Hg (25"). Pumping rate is 15 ml per stroke. Built in vacuum release valve. Can be used as pressure pump by transferring release valve cap to vacuum port. Nozzle fits standard 6 mm (1/4") i.d. tubing.

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ANALYTICAL FUNNELS

PP

Designed specifically for analytical chemistry. Each size takes a standard size filter paper. Body of funnel is exact 60° angle. Provided with external and internal ribs. Autoclavable.

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LONG-STEM FUNNELS

PP

The 60° angles sides are provided with ribs for fast filtration and allow the use of standard size filter papers.

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VISCOUS LIQUID PUMPS

PP

Dispensing pump made of PP operated by a unique lever-action stroke. Delivery of approximately 10 ounces (ab. 300 ml) per stroke, with capability of 6 gallons per minute (ab. 22-23 litres/min). Will handle the most viscous materials (equivalent to SAE 90 oil). Spigot orientation of 360° allows versatility in dispensing. Complete with shank.

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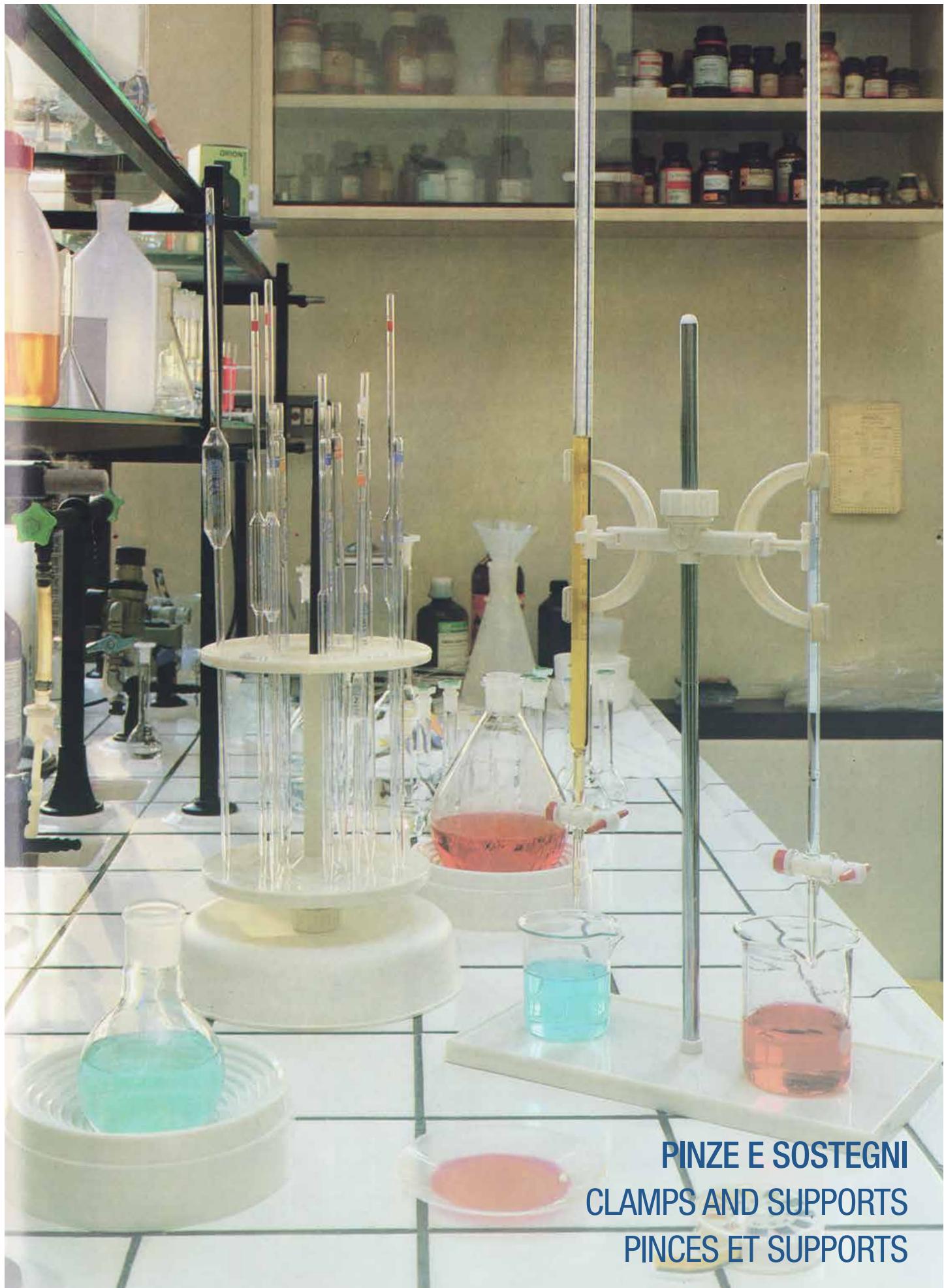
PLASTILAB

PINZE E SOSTEGNI
CLAMPS AND SUPPORTS
PINCES ET SUPPORTS

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DRAINING RACK

HIGH IMPACT PS

The back plate of the 72 place support is made from a single mould of 4 mm rigid HIPS. The peg sockets have closed ends eliminating leakage and consequent biohazards. The seamless drip channel collects waste which drains via drain tube provided. 72 push in pegs which can accommodate items with "neck" bore in excess of 15 mm are supplied and 11 small pegs of diameter 6 mm for small diameter items such as test tubes. They are easily removable for cleaning or to accommodate particularly shaped items. This rack will not stain or rust. Comes complete with mounting kit, draining hose, and fixing template. Additional racks can be joined together to make a bank of racks, using the joining plugs provided. Supplied with draining kit (connector + tubes). Discharging tube: internal 11 mm, external 13 mm.

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FUNNEL HOLDER

PP

Single or double holder for funnels as small as 25 mm (with adaptor supplied) and up to 170 mm. Without adaptor it can be used with powder funnels with stem diameter not exceeding 35 mm. Two olives provided for use with rods from 8 to 14 mm.

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SEPARATING FUNNEL HOLDER

PP

Horseshoe funnel holder especially for separating funnels. Supplied with two 'olives' so will fit supporting rod diameters 8 to 14 mm. Cannot be used with funnels below 75 mm diameter.

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CLAMPS FOR GLASS JOINTS**POM**

Plastic clamps for safe connection of standard-finish glass conical joints. Corrosion proof, do not scratch glass. Available in various colours according to dimension. Autoclavable

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GLASS FIBER SPATULAS**GLASS FIBRE FILLED NYLON**

Made according to DIN 12890. A range of robust, inexpensive general purpose spatulas. Excellent chemical resistance. Scraper plus spoon or double scraper configuration. Suitable for use in food and beverage industry. Autoclavable. Suitable for food contact. (NSF/ANSI STANDARD 51).

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Pinzette, flessibili, ad alto modulo elastico. Autoclavabili.

FORCIPS**POM**

Self-sprung forceps with rounded ends. Autoclavable.

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BURETTE CLAMPS**PP**

Boss head supplied with two olives so that the clamp can fit on to support rods having diameters 8 - 14 mm. Stainless steel spring loaded centre arm holds burette firmly against the two outer arms and is protected by rubber grips. Graduations and meniscus are visible at all times. Suitable for all burettes.

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RECTANGULAR BASE

PP WITH CHROMIUM PLATED STEEL ROD

The rods are 25 cm standard length. Rods can be screwed into one another to obtain desired height. Each base is supplied with two rods.

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ROUND BOTTOM FLASK HOLDER

PP

Round bottomed flasks from the smallest up to 10 litres can be supported safely on the internal "steps" of this support. Hot flasks up to 100°C (120°C for short periods) can be supported without damaging the holder. Autoclavable.

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BENCH PIPETTE HOLDER

PMMA

This bench pipette holder, made of particularly clear polymethylmethacrylate (PMMA), has high level finishes and its content can be seen from any angle. Complete with four tilted support shelves and white top. The four separate compartments are ideal for housing plastic or glass serologic pipettes of various capacities (from 1 ml to 50 ml). Do not shine ultraviolet light on it.

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PIPETTE STAND, VERTICAL**PP**

Accomodates 18 pipettes with diameters not exceeding 10 mm, plus 10 pipettes with diameters not exceeding 15 mm. Small holes in base allow drainage. Boss head (supplied) allows for the holder to be raised or lowered on supporting rod of 8-14 mm using the olivessupplied. Use the custom designed dished base unit with PVC rod (Art 262) to complete the system.

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**PIPETTE STAND, VERTICAL****PP**

Accomodates 8 pipettes with diameters not exceeding 10 mm, plus 8 pipettes with diameters not exceeding 15 mm. Small holes in base allow drainage.

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**PIPETTE STAND****PP**

Rotary pipettes holder 94 places. PP desktop pipette holder. Rotating base, holds up to 94 pipettes. The swivel stand accommodates up to 94 pipettes with a diameter not exceeding 14mm

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**PIPETTE JAR****PP**

Very versatile. Particularly suitable for storing pipettes. Rugged. No more chipped ends as pipettes are cushioned by PP. The jar, but not the PE cap, can be autoclaved at 120°C for 20 minutes. Tight fitting cover protects pipettes during transportation and from dust etc. Welded flanged base gives excellent stability. Can be easily cut down to suit shorter length pipettes. Can be canted on side for drainage whilst pipettes insitu.

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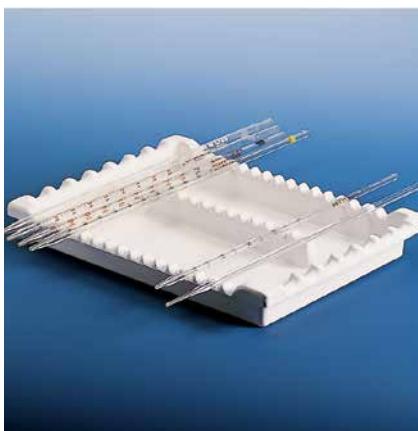
PIPETTE TRAY FOR DRAWER

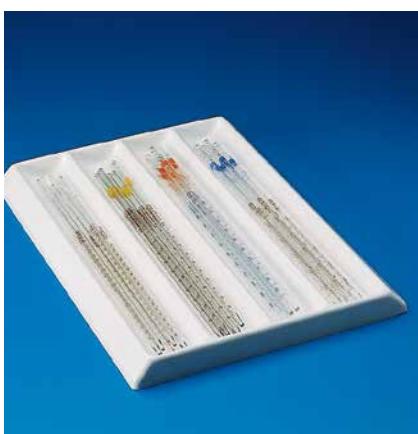
PVC

Can accommodate 7 pipettes with overall stem diameter not exceeding 20 mm longitudinally or 16 pipettes with stem diameter not exceeding 10 mm laterally. Very useful work station holder not only for pipettes.

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PIPETTE TRAY FOR DRAWER

PVC

Designed to hold about 30 pipettes. It has 4 compartments to accomodate pipettes of 1-2-5-10 ml capacities.

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PIPETTE FILLER WITH ADAPTER

RED RUBBER

A moulded rubber bulb, with three glass ball valves which can be controlled by finger pressure. There are no metal parts to corrode and the filler can be used with all liquids except those that attack rubber. Can be used on bulb or cylindrical pipettes of different capacity. Complete with adapter for pipettes with capacities more than 20 ml (up to 100 ml).

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UNIVERSAL SAFETY PIPETTE BULB

RED RUBBER AND SILICONE

Red rubber and silicone pipette bulb for pipettes up to 25ml.

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PIPETTE PUMPS

PP

Suction and dispensing is achieved by rotating the wheel. A pressure lever ensures that all liquid is evacuated.

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PIPETTE FILLER WITH PLUNGER

ABS, HDPE AND PA6

Comfortable grip. Simply press plunger and release to fill pipette. Knurled cap adjusts the plunger stroke for repeating preset volume. Rubber chuck will hold pipette securely and safely. Will handle pipettes up to 10 ml. Provided with graduations on the body. Moulded in yellow ABS.

Art.	Ø mm	h mm	Cap. max. pipette / Max. capacity pipettes / Cap. max pipettes ml
198	40	100	10

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"VADOSA" VACUUM PIPETTING DEVICE

PP

Safety vacuum pipetting device. Easy to use, being able to draw up and measure any liquid, even the most viscous ones. Vadosa fits every type and size of pipette from 0.1 ml up to 25 ml. Will deliver even the smallest quantities with simplicity, precision and speed. A simple, one handed operation with no hand squeezing, no sucking, no guessing. Made of acid and alkali-resistant PP.

Art.	Ø mm	h mm	Cap. max. pipette / Max. capacity pipettes / Cap. max pipettes ml
199	62	118	25

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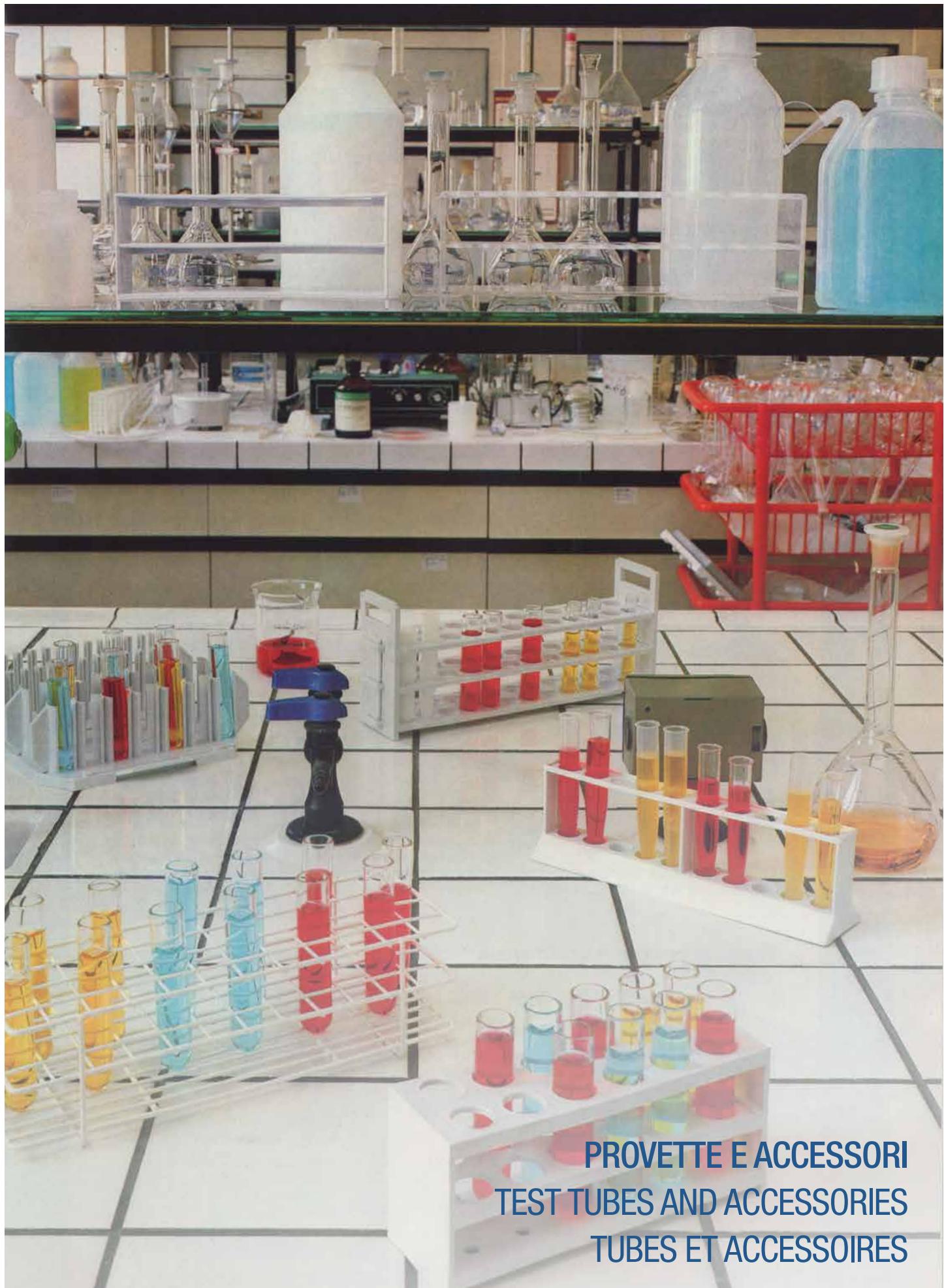
PLASTILAB

PROVETTE E ACCESSORI
TEST TUBES AND ACCESSORIES
TUBES ET ACCESSOIRES

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CENTRIFUGE TUBES,
CYLINDRICAL

PP

Autoclavable. Can withstand temperatures up to 120°C. Rimmed. High chemical resistance. Cannot be used with naked flame. Can also be used as conventional test tubes. Very good translucency.

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CONICAL CENTRIFUGE TUBES

PP

High chemical resistance; autoclavable. Can withstand temperatures up to 120°C. Cannot be used with naked flame.

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CONICAL CENTRIFUGE TUBES,
GRADUATED

PP

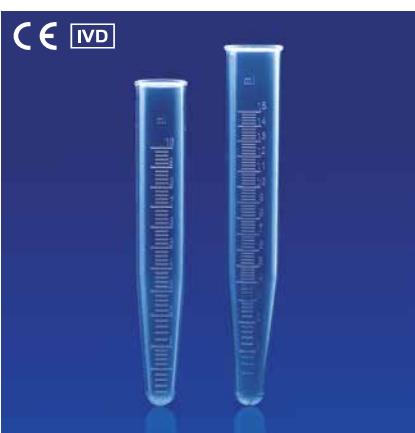
Autoclavable. Can withstand temperatures up to 120°C. Rimmed. High chemical resistance. Good transparency especially when in contact with liquids. Cannot be used with naked flame. Can be used as conventional test tubes.

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CONICAL CENTRIFUGE TUBES,
GRADUATED

PMP (TPX®)

Autoclavable. Can withstand temperatures up to 170°C. Rimmed. High chemical resistance. Crystal clear. Cannot be used with naked flame. Can be used as conventional test tubes.

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TEST TUBE CAPS

PE

Caps to fit plastic or glass test tubes.

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MICRO TEST TUBE RACK -
PREPARATION

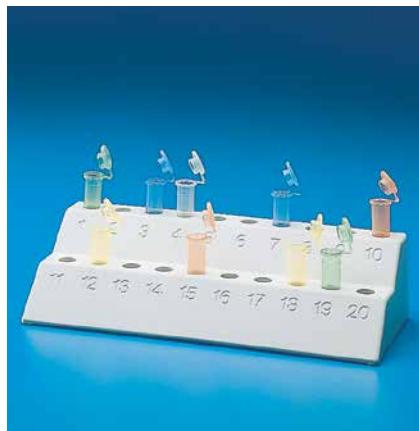
PP

Unique, dual level design for holding 20 tubes in pre-numbered cavities (1-10; 11-20). Racks have locations on base to enable attachment in rows, or back to back to each other, using optional linking clips (Art.295). Excellent for preparation work. Autoclavable.

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TEST TUBE RACKS

PP

Specifically designed for titration work. The test tubes are held in position by the upstand support. All the contents of the test tubes are visible all the time. Can withstand temperatures up to 120°C. Can be used in water baths.

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BASE PLATE FOR TEST TUBE
RACKS

PP

Base plate that can hold two of Art 130 to 135 series test tube racks in the slots provided. Useful for colorimetric analysis and to stock standard samples.

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TUBE-RACK, MICROCENTRIFUGE
(1,5 ML)

PP

Two-tier, available in four different colours. These versatile racks are autoclavable and can be used dry, in water baths or in freezers. Each rack has an alpha-numeric reference for easy sample identification. Can accommodate 100x1,5 ml tubes with closed caps or without caps, and 50x1,5 ml tubes with open caps. Racks are stackable even when loaded and have (purpose) designed label spaces on each side (for bar coding, etc.).

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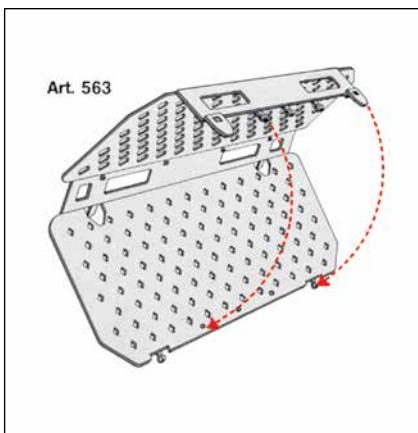
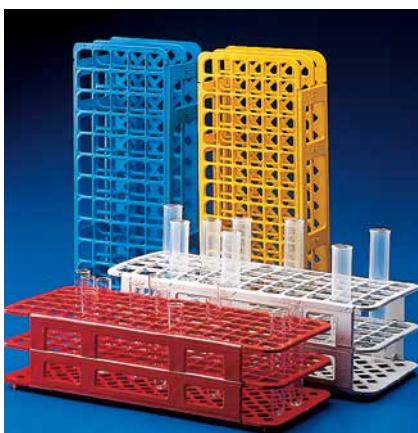
UNIVERSAL TEST TUBE RACK

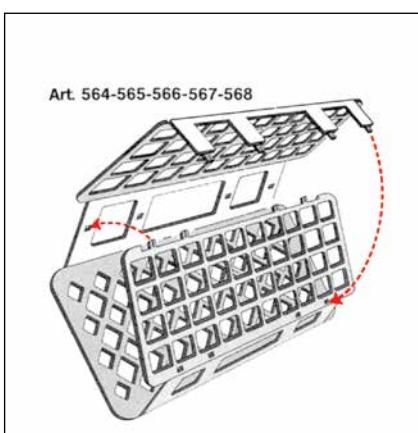
PP

Large capacity racks. Autoclavable at +121°C for 20 minutes. Can be used dry, wet, in waterbaths. Will not float. Special design allows samples to be incubated at same temperature. Can be used in freezers without deforming or going brittle. Will not rot, rust or deform. Alpha numerical grid reference on top tier for easy identification of samples. Colour coded for bath work. Stackable when empty.

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TEST TUBE RACKS TWO/THREE TIER**PP**

Lightweight, unbreakable traditional pattern test tube racks. Can withstand temperature up to 120° C. Autoclavable.

Art.	N. piani No. tiers N. étagères	Posti Places Emplac.	Ø fori mm Ø holes mm Ø trous mm	Dim. mm	h mm
559	3	12	12	190x60	80
560	3	12	20	190x60	80
562	3	24	20	375x65	95
370	2	10	18	250x60	74

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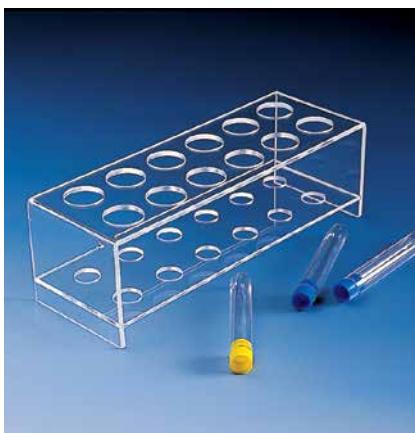
TEST TUBE RACKS TWO-TIER**PMMA (PLEXIGLAS)**

Sturdy test tube racks made of glass clear PMMA (Plexiglas). Smaller holes in the lower tier of the rack hold the test tubes safely.

Art.	Posti Places Emplac.	Ø fori super. mm Ø upper holes mm Ø supérieur trous mm	Ø fori inferiori mm Ø lower holes mm Ø inférieur trous mm	Dim. mm
889	12	32	20	90,2x255,4x81
890	12	38	22	97,5x289x81
891	12	44	22	110,5x316,5x81

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TEST TUBE TRAYS**EXPANDED PS**

Very useful trays for large quantities of test tubes. Holes are deep enough to keep test tubes, vials etc. stable. Light weight. Stackable. Finger recesses to aid lifting. 1175 will take Eppendorf® tubes type Art. 298.

Art.	Posti Places Emplac.	Ø fori mm Ø holes mm Ø trous mm	Prof. fori mm Depth of hole mm Profon. des trous mm	Dim. mm
1175	250	8,5	15	220x220x32
1177	119	14,5	18	220x220x32
1179	96	16,5	22	220x220x32

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ICE BUCKET**EXPANDED PUR**

Equipped with lid and two side-handless. Thickness of about 15 mm. Red colour. Excellent isolating feature. Can be used with dry or minced ice and cooling mixtures. Liquid nitrogen resistant for a few minutes.

Art.	Cap. lt	Ø mm	h mm
1000	4,5	270	200

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Art.
201**DRINKING STRAW****PP**

Excellent temperature resistance.
Autoclavable.

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Art.
293**TONGUE DEPRESSOR****PMMA (PLEXIGLASS)**

High mechanical resistance.
Autoclavable at 90°C.

Lungh. mm / Length mm / Long. mm
190**ABAISSE-LANGUE****PMMA (PLEXIGLASS)**

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Art.
192**INVALID CUP****PP**

Autoclavable. Removable lid for cleaning purposes.

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Art.
257**KIDNEY DISH****PP**

Excellent chemical and temperature resistance.
Autoclavable.

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Used in hospital for clearance testing on urine collected during the whole day long. Specific design with rectangular base, thought for easy storage of the bottles in the ward's cabinets and for enhanced stability when in use; the ergonomic design of the handgrip allows easy pouring. Screw capped and graduated up to 2.000 ml.

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SAMPLING BOTTLE FOR URINE

PE

Graduated up to 2500 ml. Supplied with cap for easy sampling.

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BED PAN

PP

Moulded in white PP. Can withstand temperatures up to 120°C.

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INPUT POT

PP

Smaller pot very useful in storing or wet sterilising dissecting instruments. Larger pot for longer instruments. They could also be used as drip containers or soak jars. Wide base makes them very stable. Autoclavable.

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DOUCHE CANS OR IRRIGATORS

PP

Provided with flat back and suspension tag. Autoclavable.

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FLUOROWARE

PTFE LABORATORY ARTICLES

PTFEs (Polytetrafluoroethylene) are polymers with high molecular weight that are widely used owing to their optimum chemical, electrical and thermal properties; their friction coefficient is very low. The essential features are the following.

PRACTICALLY TOTAL CHEMICAL INERTIA WITH REAGENTS AND SOLVENTS

They can be used in the more corrosive environments and do not contaminate the more sensitive materials in ultra-pure or corrosive applications.

VERY HIGH THERMAL STABILITY

Use temperature from -200°C to +300°C thermal degradation does not occur before +400°C

EXCEPTIONAL ELECTRICAL PROPERTIES

Extremely high electrical resistance low dielectric constant and low dielectric loss factor

OUTSTANDING NON-STICK PROPERTIES

Self-lubricating and shockproof features, toughness and flexibility even at low temperatures

APPLICATIONS

Chemical: very high chemical resistance to acids, alkali and solvents is typical of the fluoride polymers; it is therefore used for components in the petrochemical and chemical industry alimentary: physiologically inert, it is approved by some Authorities for uses in contact with foods electrical: excellent dielectric, self-extinguishing and stability in bad weather features make it increasingly used in this sector mechanical: its low friction coefficient makes it usable for very low load applications

The magnetic stir bars shown on the following pages have internal magnets in Alnico V that gives them of high duration and power features. These magnets are encapsulated in first quality PTFE cases. The strict quality controls applied to these products ensure correct positioning of the magnetic core, the quality of the external surface, breakage resistance and constancy of the thickness of the external covering.

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STIRRER BLADES FOR BALLOON FLASKS

PTFE

Will fit all standard button-tip stirring shafts. "Long groove" keyhole (Ø mm 9) provides easy insertion into flask neck and proper positioning for stirring action. Can be used continuously at temperatures up to 250°C and, for short intervals, to 300°C. **SUITABLE FOR FOODSTUFF.**

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WATCH GLASSES**PTFE**

Superior chemical, mechanical and heat resistance. Can be used safely for continuous service at temperatures up to 250°C. **SUITABLE FOR FOODSTUFF.**

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SUPPLE CONNECTORS**PTFE**

Specially designed to connect to misaligned glass tubes. Chemically inert. Taper conforms to DIN 12242 Standard. **SUITABLE FOR FOODSTUFF.**

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STIRRING BARS, CYLINDRICAL**MAGNET PTFE COATED**

Rounded shape, smooth walls, all-purpose. **SUITABLE FOR FOODSTUFF.**

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STIRRING BARS, MAGNETIC**MAGNET PTFE COATED**

Designed for a wide variety of applications. Savings guaranteed. Stronger mixing action at lower speeds. Can be used as an addition to any standard magnetic stir bar range, creating a viable alternative. The sizes shown below are nominal, with $\pm 5\%$ tolerance in length and $\pm 10\%$ in diameter **SUITABLE FOR FOODSTUFF.**

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STIRRING BARS, OVAL**MAGNET PTFE COATED**

Ideal for containers or bottles with round bottom.
SUITABLE FOR FOODSTUFF.

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STIRRING BALL, MAGNETIC**MAGNET PTFE COATED**

Ideal for test tubes or for eccentric stirring movements.
SUITABLE FOR FOODSTUFF.

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STIRRING BARS, CROSSHEAD**MAGNET PTFE COATED**

Special shape for small containers. Particularly effective stirring action.
SUITABLE FOR FOODSTUFF.

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STIRRING BARS, DOUBLE ENDED**MAGNET PTFE COATED**

Excellent centering, limited contact surface and high turbulence even at low speeds.
SUITABLE FOR FOODSTUFF.

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ADATTE AL CONTATTO ALIMENTARE.

STIRRING BARS, DISC**MAGNET PTFE COATED**

Convex shape to get good centring.
Particularly effective stirring action.
SUITABLE FOR FOODSTUFF.

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**TRIANGULAR STIRRING BARS,
MAGNETIC****MAGNET PTFE COATED**

Cuneiform. Ideal for dissolving solids and mixing sediments thanks to its scraping action. The angled surfaces create high turbulence even at low speeds.
SUITABLE FOR FOODSTUFF.

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ADATTE AL CONTATTO ALIMENTARE.

**STIRRING BARS, MAGNETIC WITH
PIVOT RING****MAGNET PTFE COATED**

The angled surfaces create high turbulence even at low speed. For containers with irregular bottom. The pivot ring reduces the contact surface and fosters optimal stir bar positioning.

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ADATTE AL CONTATTO ALIMENTARE.

**STIRRING BARS RETRIEVERS,
MAGNETIC****MAGNET PTFE COATED**

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LINING FOR GLASS TAPERS

PTFE

Specially designed to prevent machined glass joints from binding. Conforms to DIN 12242.

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TAPE

PTFE

Particularly useful to protect any type of screw thread. Colour: white.

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TEST TUBES

PTFE

Excellent chemical, mechanical and heat resistance. Will withstand temperatures continuously up to 250°C and for short intervals up to 300°C.

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TAPERED STOPPERS

PTFE

Excellent chemical and temperature resistance. Self-lubricating. Standard taper: 1:10. According to DIN Standard 12242.

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BEAKERS**PTFE**

Excellent chemical, mechanical and heat resistance. Will withstand temperatures continuously up to 250°C and for short intervals up to 300°C.
SUITABLE FOR FOODSTUFF.

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CRYSTALLISATION VESSELS**PTFE**

Excellent chemical, mechanical and heat resistance. Can be used safely for continuous service from -270°C up to 250°C. Particularly suitable for dissolutions and digestions with acids and corrosive agents.
SUITABLE FOR FOODSTUFF.

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PROPERTIES AND ADVANTAGES

EXCELLENT TEMPERATURE STABILITY

PFA trace analysis labware maintains its stability at temperatures from -200 °C to 260 °C (-328 °F to 500 °F).

NO LEACHABLE METALS

PFA is manufactured without metals such as calcium, aluminum, iron, magnesium, nickel, copper, manganese or zinc. These metals can leach into samples from other containers, and contaminate samples.

SUPERIOR CHEMICAL RESISTANCE

PFA is inert with most reagents, including nitric and hydrochloric acid routinely used in cleaning trace analysis labware. This stability significantly reduces the risk of cross-contamination.

EXCELLENT LONG-TERM STABILITY

PFA containers often extend the stability of low (ppb) concentration standards. Such stability reduces the time and cost of renewing trace analysis standard solutions.

HIGHLY TRANSLUCENT

Translucency is important for making accurate measurements using volumetric labware.

SIMPLIFIED CLEANING

Conventional labware for trace analysis requires time-consuming and expensive cleaning. The hydrophobic and anti-adhesive properties of the exceptionally smooth-surfaced PFA labware simplify the cleaning process, making it ideal for trace analysis.

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WIDE NECK BOTTLES

PFA

Bottles made of transparent fluorinated resin with high temperature and chemical resistance (-250°C/+270°C). Provided with leakproof closure.

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VOLUMETRIC FLASKS**PFA**

Volumetric flask made of transparent fluorinated resin with high chemical and temperature resistance (-250°C / +270°C). Supplied with standard DIN stoppers in PFA. Individually calibrated ringmark according to DIN EN ISO 1042, class "A", Standards. Laser engraved lot number.

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GRADUATED BEAKERS LOW FORM**PFA**

Transparent, graduated beakers made of PFA (Perfluoroalkoxy), a fluorinated resin with high chemical and heat resistance (-250°C/+270°C).

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SAMPLE CONTAINERS**PFA**

With screw cap (PFA) for sample collection, transport and storage.

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EVAPORATING DISHES**PFA**

With snap on lid (PE) for contamination free sample preparation.

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SAMPLE TUBE**PFA**

With calibrated ring-mark at 10 ml, and screw cap (PFA) GL 25. For sample preparation, centrifugation and for Autosampler racks (\varnothing 25 mm).

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ROUND BOTTOM FLASK**PFA**

DIN 12348 (Glass) with standard ground socket. For rotary evaporators and distillation of fine chemicals. For use under vacuum.

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DISPOSABLE CUVETTES FOR SPECTROPHOTOMETRY

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PACKING:

- From Art. 1937 to Art. 1961: 10 boxes of 100 cuvettes each (tot. 1000 cuvettes) with Kartell brand.
- From Art. 81937 to Art. 81961: 60 boxes of 100 cuvettes each (tot. 6000 cuvettes) without Kartell brand.

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CAP FOR CUVETTES

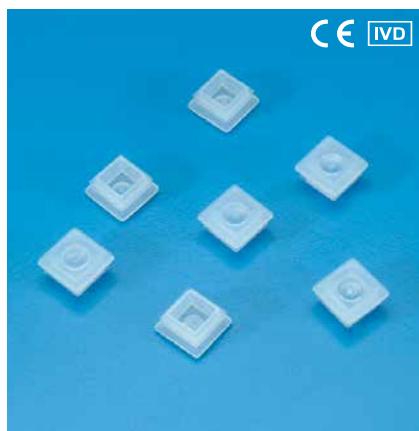
LDPE

Push-in PE stoppers for closing cuvettes with square openings (10 x 10 mm). Designed to prevent sample spillage caused by jostling and protection if sample is accidentally overturned.

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CUVETTES HOLDER

PE

Can accommodate 12x12 mm cuvettes (ext dim) cuvettes vertically in firm recesses for preparation work, or 12x10 mm cuvettes horizontally for storage and protection. In storage mode the racks are stackable.

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COBAS MIRA® MULTI-CUVETTES

PMMA

12-position polymethyl methacrylate (PMMA) multiple cuvette segments for the Cobas Mira® analyzer. A high-tech production process and stringent quality control guarantee high reproducibility. Available in racks of 15 cuvette segments for devices with continuous throughput or in single segments.

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MICRO TEST TUBES WITH CAPS FOR COBAS-BIO® AND COBAS MIRA®
PE

Micro test tubes with caps made of HDPE for Cobas Bio and Cobas Mira automatic analyzers. Available in four colours.

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MICRO TEST TUBES
PP

High transparency PP tube for easy sample checking. Flat cap with large frosted writing surface. Liquid tight for centrifugation up to RCF 21.000 x g. Large opening tab featuring rounded surfaces for smoother open/close routine and cap top with 3mm pierceable area. Autoclavable at 121°C for 20 minutes with open cap. Available in: neutral (00), blue (04), yellow (06), pink (13) and green (12).

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MICRO TEST TUBES
PP AND PE

Made of clear PP/PE they are ideal for sample analysis. It can be used for sample shipping as well. The tubes can withstand a centrifuge force of up to RCF 11.000 x g. Art. 297 and 1298 autoclavable at 121°C for 20 minutes with open cap.

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GRADUATED MICRO TEST TUBES
PP

Ideal for samples analysis. Depending on centrifuge being used, these tubes can withstand a centrifugal force of up to RCF 11.000. Flat cap with writing area. Art. 279 available in: neutral (00), blue (04), yellow (06), pink (13) and green (12).

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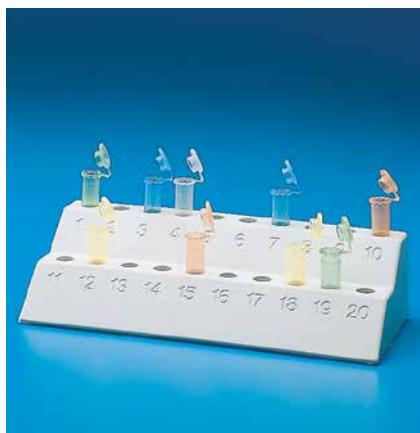
**MICRO TEST TUBE RACK
PREPARATION****PP**

Unique, dual level design for holding 20 tubes in pre-numbered cavities (1-10; 11-20). Racks have locations on the base to enable attachment in rows or back to back to each other using optional linking clips. Excellent for preparation work. Autoclavable

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Contenitori monouso in PS per contaglobuli Toa® e contaglobuli Royco-Hycel®

COULTER® CELL COUNTER CUP**PS, CAP IN PE**

Disposable PS beakers for Toa® and Royco-Hycel® cell counters

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SCINTILLATION VIALS**PE**

Unbreakable vials compatible with all liquid scintillation counters. Heavy walled to minimize solvent losses. Special PP screw cap to achieve leakproof closure (Art. 933). Insert vials, capacity of 4 ml, supplied with screw caps, (Art. 934).

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TECHNICON®, BECKMAN®, KONELAB® SAMPLE CUP
PS

PS sample cup for Technicon®, Beckman® and Konelab® analyzers. On some models an optional push-in stopper is available.

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CENTRIFICHEM®, TECHNICON®, BECKMAN® SAMPLE CUP
PS

PS sample cup for Centrifichem®, Technicon® and Beckman® analyzers. An optional push-in stopper is available for some models.

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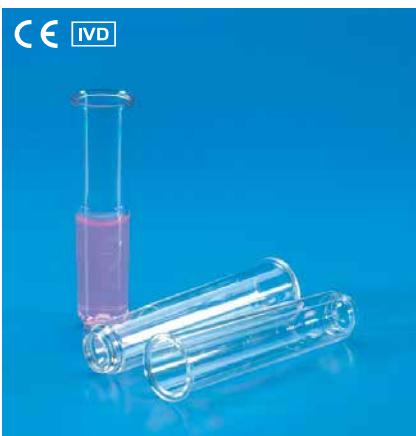
SAMPLE CUP STOPPERS
PE

Available in the anti-evaporation version for Product Code 2510 and 2511 sample cups and in the push-in version for Product Codes 2501 – 2502 – 2503 – 2510 and 2511

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CUVETTE FOR OLLI®
PS

Cuvettes for Olli-C-analyzer®, packed in PS dust proof boxes containing 192 pcs.

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CUVETTES FOR FIBRINTIMER®**PS**

Cuvettes for Fibrintimer® supplied with 500 metallic mixers.

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CUVETTES FOR AMELUNG®**PS**

Cuvettes for Amelung®, determination of coagulation time. Supplied with 1.000 metallic mixer.

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CYLINDRICAL TEST TUBES**PS**

Tight screw cap Taking into consideration the different applications, they are available sterile single packed, sterile multiple packed and not sterile.

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TEST TUBES WITH SCREW CAPS**PP**

Test tubes with quick closure screw caps (PE). Leak-proof, and suitable for centrifugation. The design of the cap avoids any contamination due to drops of liquid on the inner surface. The sterile packing (Art. 88304) is supplied with the screw cap fitted on the test tube. Available in: blue (04) - red (10)- yellow (06) - white (03).

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DISPOSABLE TUBES

PS AND PP

Disposable round or conical bottom tubes no rim, for laboratory use. Made of PS or PP. Can be used in centrifuge.

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DISPOSABLE TUBES, WITH RIM

PS AND PP

Disposable tubes with round or conical bottom, with rim, for laboratory use. Made of PS or PP. Can be used in centrifuge.

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FINNED PLUGS FOR DISPOSABLE TUBES

LDPE

Finned plugs for disposable tubes 11-13 mm or 15-17 mm in diameter. The finned plug has been carefully designed. Available colours: neutral (00), blue (04), red (10), green (12) and yellow (06).

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TAPPI PER PROVETTE MONOUSO

PE

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STOPPERS FOR DISPOSABLE TEST TUBES

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GRADUATED CONICAL TEST TUBES

PP

Supplied with PP tight screw cap. With red cap if sterile. Available also in the version with self-standing bottom. (Withstand up to -80°C).

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GRADUATED CONICAL CENTRIFUGE TUBES

PP

15 and 50ml graduated conical tubes in high-transparency PP with large whitemarking area. Withstands temperature extremes of -80°C to +121°C. Includes blue thread-design plug-seal cap with ribbed edge for safe and easy opening and closing. Conical bottom for total sample recovery. Easy-to-read gradations are pre-printed as follows: 15 ml tubes are graduated in 0.5 ml subdivisions from 2 to 15 ml 50 ml tubes are graduated in 2.5 ml subdivisions from 5 to 50 ml. Sterile tubes are irradiated by ionizing beta radiation. Designed for sample collection or transportation. Apyrogenic, endotoxines/free, autoclavable and carry the CE conformity mark pursuant to Directive 98/79/CE

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SPECIMEN CONTAINER

PS WITH PE LID

Specimen container made of PS with PE lid. Waterproof closure. Ideal for storage, collection and transportation of specimens.

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SPUTUM COLLECTION CONTAINER

PP

Autoclavabile PP sputum collection container with leakproof lid. Ideal for storage, collection and transportation of specimens.

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FAECES CONTAINER

STOPPER: PS SPOON: PP

Particularly designed for collection and delivery of faeces. Made of transparent PS with stopper of PP, incorporating a small spoon to facilitate sample handling.

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SCREW CAP CONTAINER

PS WITH PE SCREW CAP

Container with screw cap and spoon, 30 ml and 60 ml. Available sterile.

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SCREW CAP CONTAINER

PP

Cylindrical PP containers in PP material, wide mouth with screw cap lid. Available Sterile single packed and not sterile.

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URINE CUPS WITH SCREW CAP**PP**

60 ml PP urine cups with screw cap. Non-sterile cups have a blue screw cap and are available in multi-packs. Sterile cups have a red screw cap and are individually wrapped.

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URINE CUPS WITH SCREW CAP**PP**

150 and 200 ml PP urine cups with screw cap. Non sterile cups have a blue screw cap and are sold in multi-packs. Sterile cups have a red screw cap and a satinized marking area and are individually wrapped.

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CONTAINERS FOR URINE AND BIOLOGICAL SPECIMENS**PS WITH PE AND PP SCREW CAP**

60 ml PS container with PE screw cap and 200ml with PP screw cap.

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TEST TUBES WITH SCREW CAPS**PS WITH PE SCREW CAP**

12 and 25 ml PS tubes with PE screw caps. Tubes have a round bottom and a freestanding skirt. Sterile tubes are available individually wrapped. The 25-ml individual tubes have a self-adhesive label for marking.

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ERYTHROCYTE SEDIMENTATION RATE SYSTEM
PS / PP

The Sedi-Rate system allows a rapid and safe execution of the ESR test according to the Westergreen method. The system is composed of a tube in PP (12x86 mm) with label, complete with pink-cap containing 0,2 ml sodium citrate. The pipette in PS is silk-screened with a graduation scale from "0" to "180" and is provided with a fibre cap (allowing the blood to reach the 0 level) and with a little piston allowing a quick filling of the pipette thanks to the pressure exerted within the tube. Test procedure:

1. fill the test tube with whole blood up to the mark on the label (1 ml)
2. mix delicately by turning it over
3. insert the pipette until the blood reaches the 0 level
4. wait 60 minutes and read.

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10 place numbered rack for SEDI-RATE system

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TIPS, DISPOSABLE FOR AUTOMATIC MICROPPIPETTES

Manufactured from virgin cadmium-free PP. High clarity and excellent water-repellent properties designed to keep sample retention within the tip to a minimum. The precise centering of the orifice at the end of the tip guarantees reliable directionality of the sample when delivered. These tips are recommended for use with the Kartell automatic micropipettes but deliver excellent results even if used with the most other micropipette brands available on the market. Disposable Kartell tips come in three different package types to satisfy a variety of needs as economically as possible:

- bulk bag
- packed in racks
- refills for reloading racks

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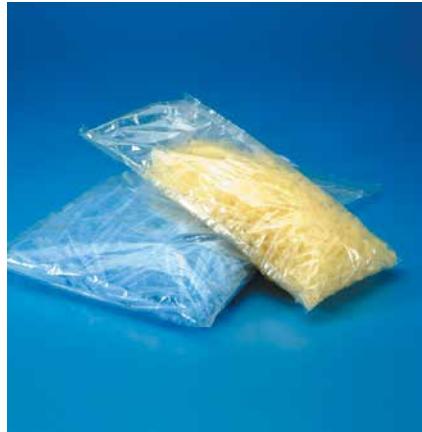
TIPS RACK

Some disposable tips are also available racked. Boxes containing the racks are in PP and have a snap-on lid. Boxes containing "macro" tips (5,000 µl and 10,000 µl) have a sliding horizontal cover. Containers and tips are autoclavable to +121°C for 20 minutes.

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TIPS REFILL

System comprising a support plate with an array of holes containing pipe tips. Designed for the reuse of empty racked tip boxes once the racked tips are used. Several tip styles are available, with the exception of the "macro" tips (5,000 µl and 10,000 µl). As an option for all refill tips, with the exception of "macro" tips (5,000 µl and 10,000 µl), spare empty racked tip boxes are available in yellow PP. As an alternative, clear polycarbonate boxes with a horizontal sliding cover can be used; there are two different formats: one for refill tips with a volume range up to 250 µl and the other for refill tips with a volume range up to 1,000 µl.

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CAPACITY 0,1-10 µL

Compatibility: Gilson®, Eppendorf®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITÉ 0,1-10 µL

Compatibilité: Gilson®, Eppendorf®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

Art.	Colore / Colour / Couleur	Tipo / Type / Type	Confezione / Packing / Conditionnement
956	Neutro / neutral / neutre	Gilson® micro	Sfusi / Bulk
9646	Neutro / neutral / neutre	Gilson® micro	Rack
9747	Neutro / neutral / neutre	Gilson® micro	Refill

**CAPACITÀ 0,5-20 µL**

Compatibilità: Kartell® Pluripet PL10, Eppendorf®, Gilson®, Biohit®, Brand®, Socorex®, Finnpipette®, Thermo®, HTL®

CAPACITY 0,5-20 µL

Compatibility: Kartell® Pluripet PL10, Eppendorf®, Gilson®, Biohit®, Brand®, Socorex®, Finnpipette®, Thermo®, HTL®

CAPACITÉ 0,5-20 µL

Compatibilité: Kartell® Pluripet PL10, Eppendorf®, Gilson®, Biohit®, Brand®, Socorex®, Finnpipette®, Thermo®, HTL®

Art.	Colore / Colour / Couleur	Tipo / Type / Type	Confezione / Packing / Conditionnement
959	Neutro / neutral / neutre	Eppendorf® Cristall	Sfusi / Bulk
9650	Neutro / neutral / neutre	Eppendorf® Cristall	Rack
9746	Neutro / neutral / neutre	Eppendorf® Cristall	Refill

**CAPACITÀ 2-20 µL**

Compatibilità: Eppendorf®, Gilson®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITY 2-20 µL

Compatibility: Eppendorf®, Gilson®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITÉ 2-20 µL

Compatibilité: Eppendorf®, Gilson®, HTL®, Nichiryo®, Finnpipette®, Thermo®

Art.	Colore / Colour / Couleur	Tipo / Type / Type	Confezione / Packing / Conditionnement
979	Neutro / neutral / neutre	Eppendorf®	Sfusi / Bulk
9657	Neutro / neutral / neutre	Eppendorf®	Rack
9745	Neutro / neutral / neutre	Eppendorf®	Refill

**CAPACITÀ 2-200 µL**

Compatibilità: Gilson®, Eppendorf®, Socorex®, Nichiryo®, Biohit®, Brand®, Hamilton Kartell® multipet

CAPACITY 2-200 µL

Compatibility: Gilson®, Eppendorf®, Socorex®, Nichiryo®, Biohit®, Brand®, Hamilton Kartell® multipet

CAPACITÉ 2-200µL

Compatibilité: Gilson®, Eppendorf®, Socorex®, Nichiryo®, Biohit®, Brand®, Hamilton Kartell® multipet

Art.	Colore / Colour / Couleur	Tipo / Type / Type	Confezione / Packing / Conditionnement
960	Giallo / Yellow / Jaune	Universal	Sfusi / Bulk
980	Giallo grad. / Yellow grad. / Jaune grad.	Universal	Sfusi / Bulk

CAPACITÀ 5-200 µL

Compatibilità: Kartell®, Eppendorf®, Gilson®, Biohit®, Brand®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITY 5-200 µL

Compatibility: Kartell®, Eppendorf®, Gilson®, Biohit®, Brand®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITÉ 5-200µL

Compatibilité: Kartell®, Eppendorf®, Gilson®, Biohit®, Brand®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

Art.	Colore / Colour / Couleur	Tipo / Type / Type	Confezione / Packing / Conditionnement
9641	Giallo / Yellow / Jaune	Universal	Rack
9749	Giallo / Yellow / Jaune	Universal	Refill

CAPACITÀ 5-200 µL**Compatibilità:** Gilson®, Kartell® multipet**CAPACITY 5-200 µL****Compatibility:** Gilson®, Kartell® multipet**CAPACITÉ 5-200 µL****Compatibilité:** Gilson®, Kartell® multipet

Art.	Colore / Colour / Couleur	Tipo / Type / Type	Confezione / Packing / Conditionnement
958	Giallo / Yellow / Jaune	Gilson®	Sfusi / Bulk

**CAPACITÀ 2-200 µL****Compatibilità:** Kartell® Multipet, Gilson®, Eppendorf®, Biohit®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®**CAPACITY 2-200 µL****Compatibility:** Kartell® Multipet, Gilson®, Eppendorf®, Biohit®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®**CAPACITÉ 2-200 µL****Compatibilité:** Kartell® Multipet, Gilson®, Eppendorf®, Biohit®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

Art.	Colore / Colour / Couleur	Tipo / Type / Type	Confezione / Packing / Conditionnement
940	Neutro grad. / neutral grad. / neutre grad.	Gilson®	Sfusi / Bulk
9643	Neutro grad. / neutral grad. / neutre grad.	Gilson®	Rack
9751	Neutro grad. / neutral grad. / neutre grad.	Gilson®	Refill

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FINE TIP 200 µL

To be used whenever it is difficult to reach the test tubes bottom with the standard products. Sterile rack of 96 fine tips 200 µl, packed conceived for Gel Loading.

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Compatibilità: Kartell®, Biohit®, Gilson®

CAPACITY 2-300 µL

Compatibility: Kartell®, Biohit®, Gilson®

CAPACITÉ 2-300 µL

Compatibilité: Kartell®, Biohit®, Gilson®

Art.
971
9644
9752Colore / Colour / Couleur
Neutro / neutral / neutre
Neutro / neutral / neutre
Neutro / neutral / neutreTipo / Type / Type
Biohit®
Biohit®
Biohit®Confezione / Packing / Conditionnement
Sfusi / Bulk
Rack
Refill

CAPACITÀ 20-300 µL

Compatibilità: Eppendorf®, Gilson®, Biohit®, Brand®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITY 20-300 µL

Compatibility: Eppendorf®, Gilson®, Biohit®, Brand®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITÉ 20-300 µL

Compatibilité: Eppendorf®, Gilson®, Biohit®, Brand®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

Art.
935
9647
9754Colore / Colour / Couleur
Neutro grad. / neutral grad. / neutre grad.
Neutro grad. / neutral grad. / neutre grad.
Neutro grad. / neutral grad. / neutre grad.Tipo / Type / Type
Eppendorf®
Eppendorf®
Eppendorf®Confezione / Packing / Conditionnement
Sfusi / Bulk
Rack
Refill

CAPACITÀ 100-1000 µL

Compatibilità: Kartell®, Gilson®, Brand®, Nichiryo®, Biohit®, Socorex®, Eppendorf®, Hamilton

CAPACITY 100-1000 µL

Compatibility: Kartell®, Gilson®, Brand®, Nichiryo®, Biohit®, Socorex®, Eppendorf®, Hamilton

CAPACITÉ 100-1000 µL

Compatibilité: Kartell®, Gilson®, Brand®, Nichiryo®, Biohit®, Socorex®, Eppendorf®, Hamilton

Art.
961Colore / Colour / Couleur
Blu / blue / bleuTipo / Type / Type
UniversalConfezione / Packing / Conditionnement
Sfusi / Bulk

CAPACITÀ 50-1000 µL

Compatibilità: Kartell®, Eppendorf®, Gilson®, Biohit®, Brand®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITY 50-1000 µL

Compatibility: Kartell®, Eppendorf®, Gilson®, Biohit®, Brand®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITÉ 50-1000 µL

Compatibilité: Kartell®, Eppendorf®, Gilson®, Biohit®, Brand®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

Art.
9640
9748Colore / Colour / Couleur
Blu / blue / bleu
Blu / blue / bleuTipo / Type / Type
Universal
UniversalConfezione / Packing / Conditionnement
Rack
Refill

CAPACITÀ 50-1000 µL

Compatibilità: Kartell®, PL1001, Gilson®, Eppendorf®, Biohit®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITY 50-1000 µL

Compatibility: Kartell®, PL1001, Gilson®, Eppendorf®, Biohit®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITÉ 50-1000 µL

Compatibilité: Kartell®, PL1001, Gilson®, Eppendorf®, Biohit®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CE IVD

Art.	Colore / Colour / Couleur	Tipo / Type / Type	Confezione / Packing / Conditionnement
941	Blu / blue / bleu	Gilson®	Sfusi / Bulk
9645	Blu / blue / bleu	Gilson®	Rack
9753	Blu / blue / bleu	Gilson®	Refill

CAPACITÀ 200-1000 µL

Compatibilità: Kartell®, Gilson®, Eppendorf®, Biohit®, Hamilton®

CAPACITY 200-1000 µL

Compatibility: Kartell®, Gilson®, Eppendorf®, Biohit®, Hamilton®

CAPACITÉ 200-1000 µL

Compatibilité: Kartell®, Gilson®, Eppendorf®, Biohit®, Hamilton®

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CAPACITY 101-1000 µL

Compatibility: Oxford®, Finnpipette®, Labsystem®

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CAPACITY 5-1000 µL

Compatibility: Beckman®, Sherwood®, Lancer®

CAPACITÉ 5-1000 µL

Compatibilité: Beckman®, Sherwood®, Lancer®

Art.
967Colore / Colour / Couleur
Blu / blue / bleuTipo / Type / Type
Beckman®Confezione / Packing / Conditionnement
Sfusi / bulk**CAPACITÀ 1000-5000 µL**

Compatibilità: Kartell®, Gilson®, Socorex®, HTL®, Nichiryo®

CAPACITY 1000-5000 µL

Compatibility: Kartell®, Gilson®, Socorex®, HTL®, Nichiryo®

CAPACITÉ 1000-5000 µL

Compatibilité: Kartell®, Gilson®, Socorex®, HTL®, Nichiryo®

Art.
984Colore / Colour / Couleur
Blu chiaro / blue light / bleuTipo / Type / Type
Kartell® Pluripet std and Pluripet II,
Gilson®, Socorex®, Nichiryo®Confezione / Packing / Conditionnement
Sfusi / bulkArt.
9648Colore / Colour / Couleur
Blu chiaro / blue light / bleuTipo / Type / Type
Kartell® Pluripet std and Pluripet II,
Gilson®, Socorex®, Nichiryo®Confezione / Packing / Conditionnement
Rack**CAPACITÀ 1000-5000 µL**

Compatibilità: Kartell®, Eppendorf®, Biohit®, Socorex®

CAPACITY 1000-5000 µL

Compatibility: Kartell®, Eppendorf®, Biohit®, Socorex®

CAPACITÉ 1000-5000 µL

Compatibilité: Kartell®, Eppendorf®, Biohit®, Socorex®

Art.
964Colore / Colour / Couleur
Neutro / neutral / neutreTipo / Type / Type
Kartell® Pluripet std, Eppendorf®,
Biohit®Confezione / Packing / Conditionnement
Sfusi / bulkArt.
9649Colore / Colour / Couleur
Neutro / neutral / neutreTipo / Type / Type
Kartell® Pluripet std, Eppendorf®,
Biohit®Confezione / Packing / Conditionnement
Rack

CAPACITÀ 500-5000 µL

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CAPACITY 500-5000 µL

Compatibility: Brand®, Finpipette®

CAPACITÉ 500-5000 µL

Compatibilité: Brand®, Finpipette®

Art.	Colore / Colour / Couleur	Tipo / Type / Type	Confezione / Packing / Conditionnement
942	Neutro / neutral / neutre	Finnpipette®	Sfusi / bulk
9847	Neutro / neutral / neutre	Finnpipette®	Rack

CAPACITÀ 2000-10000 µL

Compatibilità: Kartell®, Gilson®, Brand®, Socorex®, HTL®, Finpipette®

CAPACITY 2000-10000 µL

Compatibility: Kartell®, Gilson®, Brand®, Socorex®, HTL®, Finpipette®

CAPACITÉ 2000-10000 µL

Compatibilité: Kartell®, Gilson®, Brand®, Socorex®, HTL®, Finpipette®

Art.	Colore / Colour / Couleur	Tipo / Type / Type	Confezione / Packing / Conditionnement
943	Neutro / neutral / neutre	Gilson®	Sfusi / bulk
9848	Neutro / neutral / neutre	Gilson®	Rack

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EMPTY BOXES FOR REFILL

Empty yellow PP boxes with snap-on covers for use with refill tips with capacity up to 1000 µL.

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SEMITRSPARENT EMPTY
BOXES FOR REFILLS

Empty semitransparent boxes in polycarbonate with horizontal sliding cover. Available in short format for tips with a volume range up to 250 µL (Product Code 9755/L).

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CAPACITY 0,5-10 µL

PC/PP

Compatibility: Kartell®, Eppendorf®, Socorex®, Hamilton®. PC Rack containing PP tips.

CAPACITÉ 0,5-10 µL

PC/PP

Compatibilité: Kartell®, Eppendorf®, Socorex®, Hamilton®. Rack en PC contenant cônes en PP.

Art.	Colore Colour Couleur	Materiale rack Rack Material Matière du rack	Tipo Type Type	Confezione Packing Conditionnement
9624	Neutro / neutral / neutre	PP	Eppendorf® Cristall	Rack

CAPACITÀ 2-200 µL

PC/PP

Compatibilità: Art.9585: Gilson®, Art.9605: Kartell®, Eppendorf®, Socorex®, Finnpipette®, Brand®, Hamilton®

CAPACITY 2-200 µL

PC/PP

Compatibility: Art.9585: Gilson®, Art.9605: Kartell®, Eppendorf®, Socorex®, Finnpipette®, Brand®, Hamilton®

CAPACITÉ 2-200 µL

PC/PP

Compatibilité: Art.9585: Gilson®, Art.9605: Kartell®, Eppendorf®, Socorex®, Finnpipette®, Brand®, Hamilton®

Art.	Colore Colour Couleur	Materiale rack Rack Material Matière du rack	Tipo Type Type	Confezione Packing Conditionnement
9585	Giallo / yellow / jaune	PP	Gilson®	Rack
9605	Giallo / yellow / jaune	PP	Universal	Rack

CAPACITÀ 100-1000 µL

PC/PP

Compatibilità: Kartell®, Eppendorf®, Gilson®, Socorex®, Finnpipette®, Brand®, Hamilton®

CAPACITY 100-1000 µL

PC/PP

Compatibility: Kartell®, Eppendorf®, Gilson®, Socorex®, Finnpipette®, Brand®, Hamilton®

CAPACITÉ 100-1000 µL

PC/PP

Compatibilité: Kartell®, Eppendorf®, Gilson®, Socorex®, Finnpipette®, Brand®, Hamilton®

Art.	Colore Colour Couleur	Materiale rack Rack Material Matière du rack	Tipo Type Type	Confezione Packing Conditionnement
9615	Blu / blue / bleu	PP	Universal	Rack

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STERILE DISPOSABLE PIPETTE TIPS WITH FILTER

Made of PP and including a filter, these tips DNase, RNase and DNA free, ATP and metal free, cadmium free, apyrogenic and PCR inhibitor free. Sterilized by gamma radiation, they are autoclavable at +121°C for 20 minutes. Designed for use in molecular biology, cell cultures, bacteriology, virology, immunology and radioimmunology. Hydrophobic PE filters act as a barrier against aerosol contamination, preventing cross-contamination of samples and micropipettes. Tip orifices are sized to completely empty the tip in the event of over-aspiration. Although recommended for use in automatic Kartell micropipettes, these tips deliver optimum results when used with most micropipette brands available on the market.

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KARTELL FILTER TIPS COMPATIBILITY

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**CAPACITÀ 0,5-20 µL**

Compatibilità: Kartell® PL20, Eppendorf®, Gilson®, Biohit®, Brand®, Socorex®, Finnpipette®, Thermo®, HTL®

CAPACITY 0,5-20 µL

Compatibility: Kartell® PL20, Eppendorf®, Gilson®, Biohit®, Brand®, Socorex®, Finnpipette®, Thermo®, HTL®

CAPACITÉ 0,5-20 µL

Compatibilité: Kartell® PL20, Eppendorf®, Gilson®, Biohit®, Brand®, Socorex®, Finnpipette®, Thermo®, HTL®

Art.	Colore / Colour / Couleur	Tipo / Type / Type	Confezione / Packing / Conditionnement
9630	Neutro / neutral / neutre	Eppendorf® Cristal	Rack ster

**CAPACITÀ 0,1-10 µL**

Compatibilità: Gilson®, Eppendorf®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITY 0,1-10 µL

Compatibility: Gilson®, Eppendorf®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITÉ 0,1-10 µL

Compatibilité: Gilson®, Eppendorf®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

Art.	Colore / Colour / Couleur	Tipo / Type / Type	Confezione / Packing / Conditionnement
9638	Neutro / neutral / neutre	Gilson®	Rack ster

**CAPACITÀ 2-20 µL**

Compatibilità: Kartell® PL100, Eppendorf®, Gilson®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITY 2-20 µL

Compatibility: Kartell® PL100, Eppendorf®, Gilson®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITÉ 2-20 µL

Compatibilité: Kartell® PL100, Eppendorf®, Gilson®, HTL®, Nichiryo®, Finnpipette®, Thermo®

Art.	Colore / Colour / Couleur	Tipo / Type / Type	Confezione / Packing / Conditionnement
9639	Neutro / neutral / neutre	Eppendorf®	Rack ster

**CAPACITÀ 2-30 µL**

Compatibilità: Gilson®, Eppendorf®, Biohit®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITY 2-30 µL

Compatibility: Gilson®, Eppendorf®, Biohit®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITÉ 2-30 µL

Compatibilité: Gilson®, Eppendorf®, Biohit®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

Art.	Colore / Colour / Couleur	Tipo / Type / Type	Confezione / Packing / Conditionnement
9634	Neutro / neutral / neutre	Gilson®	Rack ster

CAPACITÀ 2-100 µL

Compatibilità: Kartell® PL100, Gilson®, Eppendorf®, Biohit®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITY 2-100 µL

Compatibility: Kartell® PL100, Gilson®, Eppendorf®, Biohit®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITÉ 2-100 µL

Compatibilité: Kartell® PL100, Gilson®, Eppendorf®, Biohit®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

Art.	Colore / Colour / Couleur	Tipo / Type / Type	Confezione / Packing / Conditionnement
9635	Neutro / neutral / neutre	Gilson®	Rack ster

**CAPACITÀ 5-100 µL**

Compatibilità: Kartell® PL100, Eppendorf®, Gilson®, Biohit®, Brand®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITY 5-100 µL

Compatibility: Kartell® PL100, Eppendorf®, Gilson®, Biohit®, Brand®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITÉ 5-100 µL

Compatibilité: Kartell® PL100, Eppendorf®, Gilson®, Biohit®, Brand®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

Art.	Colore / Colour / Couleur	Tipo / Type / Type	Confezione / Packing / Conditionnement
9632	Neutro / neutral / neutre	Eppendorf®	Rack ster

**CAPACITÀ 2-200 µL**

Compatibilità: Kartell® PL200, Gilson®, Eppendorf®, Biohit®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITY 2-200 µL

Compatibility: Kartell® PL200, Gilson®, Eppendorf®, Biohit®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITÉ 2-200 µL

Compatibilité: Kartell® PL200, Gilson®, Eppendorf®, Biohit®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

Art.	Colore / Colour / Couleur	Tipo / Type / Type	Confezione / Packing / Conditionnement
9629	Neutro / neutral / neutre	Universal	Rack ster

**CAPACITÀ 20-300 µL**

Compatibilità: Eppendorf®, Gilson®, Biohit®, Brand®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITY 20-300 µL

Compatibility: Eppendorf®, Gilson®, Biohit®, Brand®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITÉ 20-300 µL

Compatibilité: Eppendorf®, Gilson®, Biohit®, Brand®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

Art.	Colore / Colour / Couleur	Tipo / Type / Type	Confezione / Packing / Conditionnement
9637	Neutro / neutral / neutre	Eppendorf®	Rack ster

**CAPACITÀ 50-1000 µL**

Compatibilità: Kartell®, PL1001, Gilson®, Eppendorf®, Biohit®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITY 50-1000 µL

Compatibility: Kartell®, PL1001, Gilson®, Eppendorf®, Biohit®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITÉ 50-1000 µL

Compatibilité: Kartell®, PL1001, Gilson®, Eppendorf®, Biohit®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

Art.	Colore / Colour / Couleur	Tipo / Type / Type	Confezione / Packing / Conditionnement
9633	Neutro / neutral / neutre	Eppendorf®, Gilson®	Rack ster.

**CAPACITÀ 50-1000 µL**

Compatibilità: Eppendorf®, Gilson®, Biohit®, Brand®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITY 50-1000 µL

Compatibility: Eppendorf®, Gilson®, Biohit®, Brand®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

CAPACITÉ 50-1000 µL

Compatibilité: Eppendorf®, Gilson®, Biohit®, Brand®, Socorex®, HTL®, Nichiryo®, Finnpipette®, Thermo®

Art.	Colore / Colour / Couleur	Tipo / Type / Type	Confezione / Packing / Conditionnement
9631	Neutro / neutral / neutre	Eppendorf®	Rack ster.

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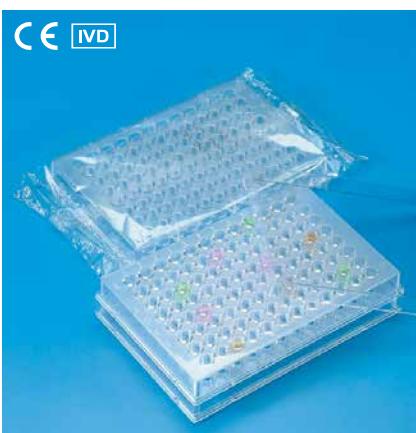
EMPTY BOXES FOR RACKED PIPETTE FILTER TIPS

Gray empty PP boxes with snapon lid for pipette filter tips.

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Multiple-well plates with 96 sample wells in clear PS. Available in three shapes: flat bottom, U bottom and V bottom wells. Aseptic and single-wrapped. PS covers are sold separately. Available also in sterile version.

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PASTEUR PIPETTE

PE

Disposable PE pipette with bulb and graduations. The slender tip guarantees precision transfer of the sample. Non-sterile pipettes are available in multi-packs; sterile pipettes are individually wrapped.

Art.	Lungh. TOT. mm TOT. Length mm Long. TOT. mm	Ø Stelo mm Stem Ø mm Tube Ø mm	Gocce/ml Drops/ml Gouttes/ml	Vol. Bulbo Bulb Draw Vol. Bulbe	Grad. ml	Grad. intvl.	Conf. Pz. Pkg.pcs Pqt.pcs
329	150	2,5	28	3,5 ml	-	-	Sing. Steril.
330	150	5,0	23	3,5 ml	1	0,25 ml	Sing. Steril.
334	150	7,8	21	3,5 ml	3	0,5 ml	Sing. Steril.
88335	150	2,5	28	3,5 ml	-	-	500
88336	150	5,0	23	3,5 ml	1	0,25 ml	500
88337	150	7,8	21	3,5 ml	3	0,5 ml	500

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DISPOSABLE DROPPING PIPETTES - BELLOW TYPE

PE

Blow-moulded in one piece with high capacity bellows reservoir for take up and expulsion. Ideal replacement of the traditional glass Pasteur pipette and rubber teat. They are graduated to give an indication of the amount of liquid being taken up and have a strong suction capacity due to the bellow-shaped leakproof bulb.

Art.	Lungh. mm Length mm Long. mm	Cap. ml	Ø stelo mm Stem Ø mm Tube Ø mm	Grad. ml
315	134	1,5	8	0,5
316	195	5,0	10	1,0

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RODS FOR CLOT DETACHMENT AND EXTRACTION

HIGH IMPACT PS

Disposable shock-resistant PS rods for detaching and extracting clots.

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INOCULATING LOOPS AND NEEDLES

HIGH IMPACT PS

Sterile inoculating loops in polybags of 20 pcs. Neutral type has a 1 µl loop at one end and a needle at the other. Blue type has a 10 µl loop at one end and a needle at the other.

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L-SHAPED SPREADERS**PS**

Disposable L-shaped sterile PS spreaders for bacteriology with special 100° angling for easier use. Especially handy for dispersing liquids across agar culture plates. Available single-wrapped or in packs of 5.

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DISPOSABLE PETRI DISHES

Disposable planar, clear Petri dishes in PS. Available in several diameters, ventilated and unventilated, aseptic or radiation- sterilized. Styles with 2, 3 and 4 separate compartments. Designed for routine lab work, bacteriological research. CE Marked.

- With aeration notches on the lid to prevent condensation formation.
- Without aeration notches to avoid agar desiccation after incubation. Gas exchange outside Petri dish is reduced thanks to the absence of notches.

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METAL TRAYS FOR HISTOLOGY**STAINLESS STEEL**

Made of stainless steel for easy removal of paraffin inclusions: they allow a precise location of the specimen into the paraffin. Packed in bags of 10 pieces.

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Art.
2922-04

TISSUE EMBEDDING SPONGE**SYNTHETIC SPONGE**

Synthetic sponge with a calibrated porosity that can easily be inserted into the tissue embedding cassettes. Very useful for small specimens in biopsy tests.

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Art.
2925

TISSUE EMBEDDING CASSETTES MEGA TYPE**POM**

Tissue embedding cassettes Mega type for medium to large-sized samples. Available in the colours: White (03), Blue (04), Yellow (06), Green (12), Pink (13).

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Art.
2919

EMBEDDING RINGS**ABS**

Embedding rings made from ABS. Available in the colours: White (03), Blue (04), Yellow (06), Green (12), Pink (13).

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Art.
2921

TISSUE EMBEDDING CASSETES, UNI TYPE**POM**

Tissue embedding cassettes for paraffin inclusions, Universal type, provided with clipon closure lid that can easily be removed. Available in the colours: White (03), Blue (04), Yellow (06), Green (12), Pink (13), Red (10).

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BALANCE BOATS

HIGH IMPACT PS

Diamond (DS) and square shaped (SQ) lightweight weighing boats provide an easy and inexpensive system for weighing any medium. They offer the following benefits:

- non contaminating (made of HIPS with hydrophobic surface)
- easily fashioned to make a funnel shape
- the DS model is available in either white (03) or black (09) so the operator can easily see that no trace of sample remains adhering to the boat
- reproducible tare weight: batch to batch variation is normally less than 5%
- **Antistatic**

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**Technical specifications common to the three series:****DOSIPET**

Dosipet is a series of adjustable volume repetitive dosing devices. They are made of high quality materials ensuring considerable mechanical and chemical resistance. They are available in three different versions, depending on the reagents to be dispensed. Possibility to choose between:

DOSIPET B

For universal use; ideal for use with reagent of average aggressiveness.

DOSIPET S

Useful for dosing concentrated acids, polar and no-polar solvents.

DOSIPET HF

Especially developed for hydrofluoric acid dosing. They are supplied complete with tubes for filling/discharging, as well as with a set of adaptors: for Dosipet 2,5/5 and 10 ml adaptors A25, A28, A45, S40; for Dosipet 25/50 and 100 ml adaptors A28, A32/45, A40.

Technical specifications common to the three series:

- precise and quick volume adjustment using the analogue slider
- precise graduated scale to ensure reproducibility of the dispensed volumes
- no loss of reagent when eliminating air or filling
- dispensing without bubbles forming
- the piston completely transfers the liquid with every movement
- total recovery of the liquid in the discharge tube by turning the delivery tip 180°
- autoclavable at +121°C for 20 minutes
- EC conformity

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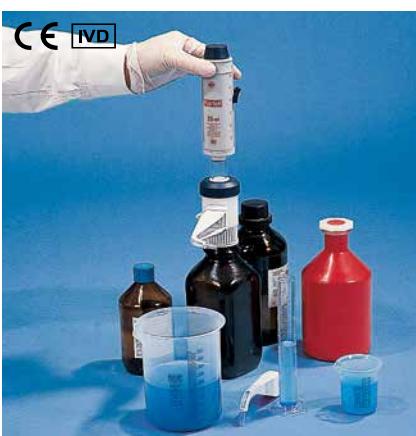
DOSIPET B

Series that can be used with every water solution type (except those containing hydrofluoric acid), marked by a high degree of convenience and reliability. It is equipped with a 3.3 borosilicate glass cylinder and PTFE piston

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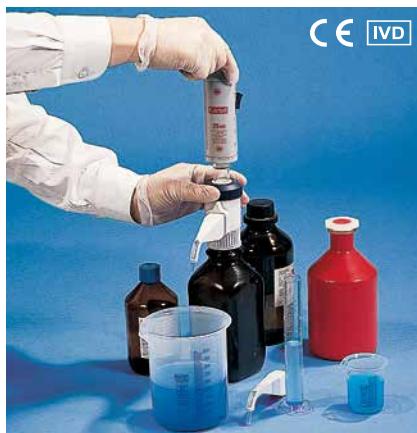
DOSIPET S

High quality components and assembly make this series the ideal choice for measuring out organic polar and non-polar solvents (see technical manual), including halogenated solvents, and concentrated mineral acids, including particularly volatile acids such as nitric and trifluoropropiatic acids. In addition, they can be used with peroxides, hydrochloric acid and sulphuric acid. It comes with a borosilicate glass piston. The reagents come into contact only with parts made of PTFE.

Art.	Mod.	Cap. ml	Subdiv. ml	Toll. ml	V ml	Filettatura GL / Thread GL / Filetage GL
83070	Dosipet S-10	1,0 - 10,0	0,20	± 0,06	0,010	32
83071	Dosipet S-25	2,5 - 25,0	0,50	± 0,15	0,025	45
83072	Dosipet S-50	5,0 - 50,0	1,00	± 0,30	0,050	45

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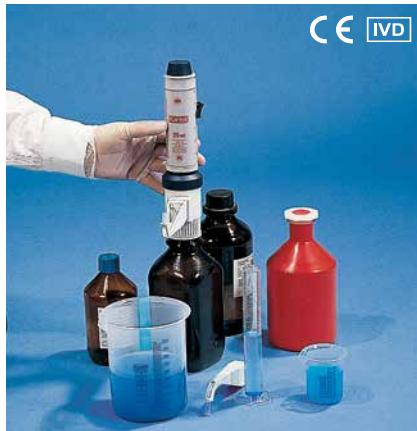
DOSIPET HF

These dosing devices are designed for specific use with hydrofluoric acid solutions. The internal components that come into contact with these solutions are made with high quality and high resistance materials (PTFE and platinum-iridium), completely inert with hydrofluoric acid. Complete with PTFE piston

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DOSILAB

UPGRADE - DosiLab is a series of adjustable volume bottle top dispenser for dosing chemicals with precision, accuracy and inexpensiveness. It is totally safe for operators and environment.

The six models of the line (0,25 ml to 100 ml) are simple to use and easy to disassemble to perform cleaning and maintenance operations; autoclavable without need for disassembly.

All the internal components that during use are in contact with liquid solutions are made of PTFE, FEP and borosilicate glass. They therefore possess excellent compatibility with all reagents, except for hydrofluoric acid. Volume adjustment is made precise and quick by using the adjustment knob, furtherly improved and supplied with internal and external locking system, to avoid accidental displacement during use. Furthermore, the dispenser allows a 360 ° rotation when placed on the bottle. A careful choice of raw materials, as well as several quality controls carried out during production process guarantee DosiLab's high precision and accuracy. Every single appliance is individually calibrated according to the ISO 8655 standards and it is supplied with its own calibration certificate.

The supply includes five adaptors for use with the most common reagent bottles (28, 32, 36, 40 and 45 mm) and a practical device for easy recalibrations that can be directly performed in laboratory, in conformity with GLO/ISO standards

TECHNICAL SPECIFICATIONS

- Glass fiber polypropylene body to increase the dimensional accuracy and the robustness of the dispenser
- Completely autoclavable at +121°C without need for disassembly.
- PTFE piston with silicone O-ring ensuring a high degree of accuracy.
- Internal cylinder in borosilicate glass protected by PP sheath.
- Valve housing: PTFE body in a single piece ensuring high chemical resistance.
- Adjustable dispensing spout with closing cap to prevent accidental dripping and contact with dangerous chemical substances
- Accuracy & Reproducibility: according to the ISO 8655 standards
- Compatibility: excellent with all reagents including organic solvents and concentrated acids, except hydrofluoric acid (HF).
- Recalibration: simple and quick with its special accessory.

Comparison of the specifications required by the ISO 8655 standards and the DOSILab specifications

The following table certifies that the DosiLab specifications fully meet the requirements required by ISO 8655, international standard for piston-driven volumetric appliances.

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Alkali	Inorganic Solutions	Acids	Organic Solvents		Hydrofluoric Acid (HF)
			Polar	Non-polar	
DosiLab					

Art.	Mod.	Cap. ml	Subdiv. ml	Prec.ml	CV ml	Flettatura GL / Thread GL / Filetage GL
83082	DosiLab 2	0,2 - 2,5	0,05	0,015	0,005	30
83083	DosiLab 5	0,5 - 5	0,10	0,025	0,005	30
83084	DosiLab 10	1 - 10	0,20	0,050	0,010	30
83085	DosiLab 30	2,5 - 30	0,50	0,150	0,030	30
83086	DosiLab 60	5 - 60	1,00	0,300	0,060	30
83087	DosiLab 100	10 - 100	2,00	0,500	0,100	30

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TECHNOTRATE DIGITAL BURETTE

Technotrate is a digital burette ideal for exact titrations directly from bottles, both in the laboratory and in the field. Developed in compliance with EN ISO 9001, it is EC certified. It has a plastic body and a large liquid crystal display also for continuous display of the dispensed volume. It adequately allows dosage volumes of up to 50 ml with 0.01 ml increases to be covered and the dosage velocity can be selected from fast mode to drop-by-drop mode.

Simple calibration

Permanent calibration refers to water at +20°C. In any case, when using liquids with a different viscosity/density, Technotrate can easily and quickly be recalibrated by pressing the dedicated button.

Practical and quick preselection of the titration volumes

Titration volume preselection is made particularly simple by pressing only the dedicated "DOS" button, looking at the display, and with just a few turns of the handle the burette cylinder is filled. The special memory button ("MEM") helps you avoid complex calculations.

Easy and safe

By turning 180°, the delivery tip makes all the liquid in the discharge tube flows inside the bottle and prevents the spilling of inappropriate liquids. The components that come into contact with the reagents are made of high resistance plastic.

Technotrate is tested one by one at the end of production, has quality certification and an individual serial number.

It is supplied in series with A25, A28, A32/45 and S45 threaded adaptors and PTFE aspiration, discharge and recirculation tubes

TECHNICAL SPECIFICATIONS:

- auto-off function after 3 minutes of non-use
- use temperature from +10°C to +40°C
- transport and storage no warmer than +60°C
- rechargeable battery: the intermittent indicator inside the display informs of the need for recharge without jeopardising the titration in progress
- precision mechanism for simple filling and elimination of air without reagent loss problems
- dispensing without bubbles forming
- battery charger: 100-240V/ 50-60Hz

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TECHNICAL SPECIFICATIONS

- the handle connection ensures a more comfortable and non-fatiguing grip
- the volume can be selected by turning the emptying button or turning the cylinder supplied for the purpose
- the large display ensures simple reading reading of the set volume value
- the button of each micropipette has a specific colour so identification of the usable volume interval is simple
- the tip ejector is adjustable in height to house a large variety of tips and is can be disassembled to simplify cleaning operations
- the recalibration device ensures quick and precise recalibrations
- the robust structure of the PVDF body ensures high mechanical and chemical resistances
- totally autoclavable at +121°C for 20 minutes
- the materials used for the structure are resistant and durable, and permit exposure to U.V. rays
- in accordance with EN IS 8655

MICROPIPETTE AUTOCLAVABILI A VOLUME VARIABILE - PLURIPET

Eight variable volume single-channel micropipettes make up the PLURIPET line equipped with an accurate and precise pipetting mechanism. They cover the range of capacities from 0.2 µl to 10,000 µl. The PL2 and PL10 models are ideal for molecular biology techniques. The PL20, PL100, PL200 and PL1.000 models are useful for dispensing small volumes of aqueous fluids of moderate density and viscosity. The PL5.000 and PL10.000 macropipettes are ideal for dispensing large volumes of rather thick and viscous aqueous fluids. They are equipped with a special filter that protects the internal piston by increasing its lifetime and preventing contaminations. They are supplied complete with a recalibration device and adhesive support for attachment to the laboratory bench. Individually calibrated with their certificates included in the package; EC/IVD-compliant.

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FIXED VOLUME AUTOCLAVABLE MICROPIPETTES - FIXOPET

The FIXOPET line is formed by eleven single channel micropipettes with fixed volume of high precision. The capacities available range from 5 µl to 1,000 µl. Each of them is equipped with adjustable tip ejector in order to eliminate any risk of contamination. The structural features of these micropipettes considerably reduce the force necessary for pipetting. Ideal for clinical and educational uses, compliant with EC/IVD.

TECHNICAL SPECIFICATIONS

- The special handle connection ensures a more comfortable and non-fatiguing grip
- the rotary device for the tip ejection does not create strain on the tendons of the thumb
- the removable ejector allows it to be used with test tubes having a particularly narrow opening
- the integrated recalibration system ensures quick and precise recalibrations
- the robust structure of the body made of PVDF ensures high mechanical and chemical resistances
- the materials used for the structure are resistant and durable, and permit exposure to U.V. rays
- totally autoclavable at +121°C for 20 minutes
- in accordance with EN IS 8655

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SPARE PARTS SET AND MICROCHANNEL MICROPIPETTES STAND

Spare part set:

- 1 cone
- 2 PTFE seal
- 2 O-ring
- 1 cone nut

Art.	Micropipette mod.
1356	FX 750 - FX 1000
1340	PL 2
1341	PL 10
1342	PL 20
1343	PL 100
1344	PL 200

Bench holder in transparent plastic material for 3 up to 6 Kartell micropipettes (Pluripet, Pluripet II, Fixopet, Labofix) or other brands. Strong structure, easy cleaning, wide base allowing great stability. Radiation sterilizable.

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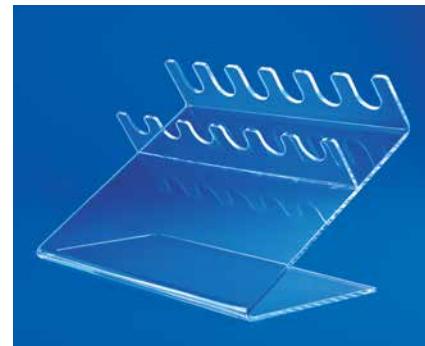
MULTICHANNEL VARIABLE VOLUME MICROPIPETTES - MULTIPET

Multichannel micropipettes with air cushion ideal for dispensing in multiwell plates for ELISA dosages, molecular biology and DNA amplifications. The special PP/PVDF structure makes them lightweight and comfortable. The pipetting mechanism makes the work very precise and the adjustment, which can be carried out with just one hand, is simple and progressive. The exclusive operation system reduces the force necessary for executing the pipetting and suction; it considerably reduces the force necessary for inserting and ejecting tips.

TECHNICAL SPECIFICATIONS

- The volume can be easily selected by turning the emptying button or turning the cylinder supplied for the purpose
- the large display ensures easy reading of the set volume value
- the MULTIPET cones are independent, so the required force for securing the 8 or 12 tips is uniformly distributed: minimum pressure suffices for inserting and ejecting them. In addition, the movement of the cones allows the micropipette to oscillate during the loading operations
- the lower part can freely turn 360°, the ejector is easily adjustable to house a large variety of tips and can be removed for cleaning
- they are totally autoclavable at +121°C for 20 minutes, without having to recalibrate them
- the materials used for the structure are resistant and durable, and permit exposure to U.V. rays
- individually calibrated with their certificates included in the package; EC/IVD-compliant
- in accordance with EN IS 8655

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FIXED VOLUME AUTOCLAVABLE MICROPIPETTES – LABOFIX

This line of micropipettes is designed to meet the increased demand for fixed volume micropipettes that combine a good quality level with a certain interesting price. Available ranging from the volume of 5 µl up to the volume of 1.000 µl. Ideal for applications in Clinical Diagnostics, Molecular Biology, Cellular Biology and Analytical Chemistry and in all those fields in which a fixed dosage carried out with good precision and accuracy is necessary. The particularly lightweight body makes the handle comfortable. Individually calibrated according to the ISO8655 standards, the new international standards for piston-driven volumetric appliances; EC/IVD-compliant.

TECHNICAL SPECIFICATIONS

- made of high quality plastic
- equipped with tip ejector and stainless steel piston for use prolonged over time
- the handle connection ensures a more comfortable and non-fatiguing grip
- the tip holder cone can house a wide variety of tips; it has a fine shape to make getting inside test tubes or containers having particularly narrow openings or lengthened shapes easier
- the robust structure of the body ensures high mechanical and chemical resistances, making them totally autoclavable at +121°C for 20 minutes

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VARIABLE VOLUME AUTOCLAVABLE MICROPIPETTES - PLURIPET II

Line made up of eight models having air displacement and equipped with coaxial ejector. Useful for applications in Molecular Biology, Microbiology, Immunology, Analytical Chemistry, etc. They ensure a high degree of precision and accuracy combined with a lower cost. The particularly light and balanced body makes the handle comfortable. They cover the range of capacities from 0.2 µl to 10,000 µl. The Pluripet II 5.000 and Pluripet II 10.000 macropipettes are ideal for dispensing large volumes of rather thick and viscous aqueous fluids. They are supplied complete with adhesive support for attachment to the laboratory bench. Individually calibrated according to the ISO8655 standards, the new international standards for piston-driven volumetric appliances; EC/IVD-compliant.

TECHNICAL SPECIFICATIONS

- made of high quality plastic and equipped with steel piston for prolonged use over time
- the handle connection ensures a more comfortable and non-fatiguing grip
- the desired volume can be set by turning the emptying button; it proves easy and safe thanks to the snap selection
- the large display ensures simple reading reading of the set volume value
- the button of each micropipetta has a specific colour so identification of the usable volume interval is simple
- the tip holder cone can house a wide variety of tips
- the recalibration device ensures quick and precise recalibrations
- the robust structure of the body ensures high mechanical and chemical resistances, making them totally autoclavable at +121°C for 20 minutes

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MINIFIX

An innovative product line including ten fixed-volume air displacement mini-micropipettes. Fully autoclavable, they are designed for fast, repetitive and precise operation and are competitively priced. Very compact in terms of size and weight, each mini-micropipette is 13 cm long and weighs approximately 10 grams. Designed for routine or biological hazard work, in situ analyses and sampling, diagnostic kits and instruction. Kartell tips fit on the conical shaft as well as most other commercially available tips. The plunger button has a different color based on dosing capacity. Individually calibrated in compliance with ISO8655, the new international standard for piston-operated volumetric apparatus. Carries the CE conformity mark.

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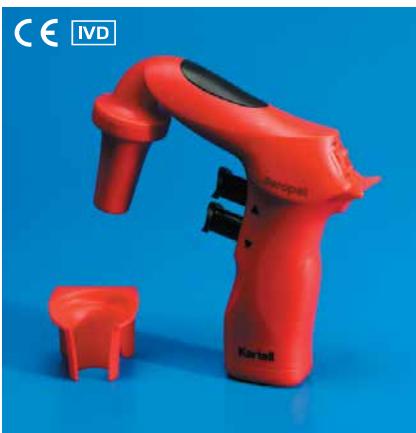
SEROPET

Seropet was developed for aspirating liquids by glass or plastic pipettes ranging from the volume of 1 ml up to 100 ml. The particularly light structure ensures comfortable use even when used for a long time and its particular configuration also makes it ideal for underhood uses. The buttons are very easy to access for choosing different modes depending on the volume pipette being used and viscosity of the liquid to aspirate. It is protected from crossed contamination in the event of excessive aspiration by the filter on the aspiration device and the safety valve. It is supplied complete with battery charger, protective PTFE hydrophobic filter (0.2 µm) and an adhesive support for attachment to the laboratory bench. EC/IVD-compliant.

TECHNICAL SPECIFICATIONS

- two different work modes: quick and normal, with continuous or pulsating adjustment
- the pipette housing is autoclavable
- supplied with safety valve to block aspiration excesses
- equipped with autoclavable protective hydrophobic anti-aerosol filter
- the special parking fins make it possible to set it down on the bench to prevent contamination of the pipette being used
- highly sensitive internal valves to get precision also with small volumes
- the NiMH battery supplied has an autonomy of up to 8 hours with continuous work and up to 1,000 charge cycles
- the indicator light warns when the battery charge is about to become exhausted

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LABOFILL - CORDLESS ELECTRONIC PIPETTE FILLING DEVICE LABOFILL

A new pipette filling device was developed to complete Kartell entry level line of liquid handling products. This device is used for filling and dispensing liquid in glass or plastic pipettes, both graduated and volumetric, up to 100 ml. LaboFill is a safe, efficient and economical instrument for liquid filling and dispensing. Due to its design is also comfortable and easy to use. LaboFill is CE marked, is individually tested after production process and is calibrated according to ISO 8655 standards.

PRODUCT FEATURES:

- Red and green L.E.D.s: the red light indicates low battery charge; green L.E.D. indicates charging status.
- Thumb wheel is provided at the back of the instrument to increase or reduce pump speed according to user's needs.
- High aspirating and flow-out speed: 25 ml in only 5 seconds.
- Silicon universal pipette housing can accommodate standard pipettes from 1 ml to 100 ml. Inside collet knurlings facilitate tight fit even with wet pipettes.
- Comfortable handgrip ensures optimal grip and prevent user from tiredness
- 0.2um PTFE hydrophobic filter is provided in the collet assembly to avoid aerosol formation and liquid entering the instrument body.
- Rechargeable battery Ni-MH 3.6 provided with the instrument, ensures cordless work for up to eight hours continuing after recharging.
- Completely battery recharging in 5 hours.
- Adaptor compatibility with 220 V and 110 V

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**ABBREVIATIONS, TEMPERATURE
AND CHEMICAL RESISTANCE OF
PLASTICS**

The table below lists commonly used abbreviations for plastics. This list covers plastics commonly employed in the manufacture of plastic laboratory ware.

Temperatures appearing in parentheses: limits tolerated for intervals only.

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**"CLEAN" AND ECOLOGICAL
PLASTICS**

Thanks to technological innovation and progressive adaptation to the most elevated eco-compatibility standards, Kartell plastic raw material are of excellent quality. In fact the used plastic materials grant wide compatibilities; hereunder you can find some standards conformities related to their characteristics.

- Foodstuff contact suitable (National and European Directive CE 1935/2004)
- Foodstuff contact suitable (US FDA CFR 170/199)
- PTFE: foodstuff contact suitable (FDA CFR TITLE 177.1550)
- Absence or limitation according the International Directives of: heavy metals, biphényles and phthalates
- RoHS Directives Conformity
- BSE and/or TSE Directives Conformity
- Atex Directives Conformity

Kartell Labware promotes the research in the field of the plastic materials, through the direct contact with the most important worldwide raw materials manufacturers and the research of innovative materials. We would like to underline that Kartell Labware materials, if not chemically contaminated, are totally recyclable

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CHEMICAL RESISTANCE

Data for the chemical resistance of salts also apply to their aqueous solutions.

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STERILISING PLASTIC LABORATORY WARE

Before sterilising any items of plastic laboratory ware, verify that no contamination or residues are present. Their presence could destroy plastics during sterilisation or autoclaving.

Observe the tolerated temperature ranges of plastic when autoclaving plastic laboratory ware. Remove any stoppers, fittings, or caps from plastic laboratory ware prior to autoclaving. Plastic vessels should be autoclaved separately from their closures and other fitting. Autoclaving vessels with their closures in place can lead to deformation and destruction of the vessels. All statements are advisory only, and imply no liability on the part of Kartell.

All statements relating to the resistances of plastic laboratory ware to high temperatures, chemicals, and to sterilisation and cleaning procedures have been cautiously formulated, based on statements of raw material manufacturers, on statements appearing in the literature, plus experience gained in actual practice.

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CLEANING PLASTIC LABORATORY WARE

All polyolefins, such as LDPE, HDPE, PP and PMP (Tpx®), as well as the fluorinated hydrocarbons PTFE, PFA, FEP, ETFE and E-CTFE, have wettable surfaces that are both highly resistant to high temperatures and chemical attack and easy to clean. Slight contamination can be removed using a chemically neutral (pH 7) cleaning agent. Heavy contamination can be removed using an alkaline (pH up to 12) cleaning agent.

Use only chemically neutral (pH7) cleaning agents on polycarbonate (PC) or Polystyrene (PS).

Laboratory ware used in trace analyses should be cleaned in a 1-N hydrochloric acid (HCl) solution for periods of not more than 6 hours and then rinsed in distilled water in order to preclude contamination by cations or anions.

Never use scouring powders or abrasive sponges when cleaning plastic laboratory ware. Use no alkaline cleaning agents on polycarbonate (PC) laboratory ware.

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A seguito di eventuali aggiornamenti, Kartell S.p.A.
After eventual revisions, Kartell S.p.A.
A la suite d'éventuelles mises à jour, Kartell S.p.A.
Kartell S.p.A.